



OceanOptics




OCEAN FLAMENIR+ SERIES

Compact Spectrometers

Installation and Operation Manual

Locations

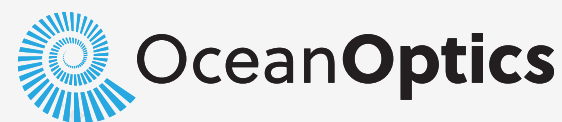

Rochester,
NY, USA

Orlando,
FL, USA

Duiven,
NL

Stuttgart
(Ostfildern), DE

Shanghai,
PRC

Americas

Manufacturing & Logistics
3500 Quadrangle Blvd.,
Orlando, FL 32817, USA

Sales: info@oceanoptics.com
Orders: orders@oceanoptics.com
Support: support@oceanoptics.com

Phone: +1 727-733-2447
Fax: +1 727-733-3962

Photonic Systems
176 Anderson Avenue, Suite F304,
Rochester, NY 14607 USA

www.oceanoptics.com

Europe, Middle East & Africa

Sales & Support
Geograaf 24, 6921 EW Duiven,
The Netherlands

Logistics
Maybachstrasse 11,
73760 Ostfildern,
Germany

Email: emeasales@oceanoptics.com

Netherlands: +31 26-319-0500
Netherlands Fax: +31 26-319-0505
Germany: +49 711-341696-0
UK: +44 1865-819-922
France: +33 442-386-588

Asia

Ocean Optics China
3F, Building 16, No. 155
Yuanke Road, Minhang District
Shanghai 200336
People's Republic of China

Email: asiамkt@oceanoptics.com

China: +86 400-623-2690
China Fax: +86 21-6295-6708
Japan & Korea: +82 10-8514-3797
Ocean Optics India: +91 81 7847-0108
Prestige Shantiniketan, Gate No. 2,
Tower C, 7th Floor, ITPL Main Road,
Whitefield, Bengaluru, Karnataka 560 048, India

Copyright © 2026 Ocean Optics.

All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted, by any means, electronic, mechanical, photocopying, recording, or otherwise, without written permission from Ocean Optics.

This manual is sold as part of an order and subject to the condition that it shall not, by way of trade or otherwise, be lent, re-sold, hired out or otherwise circulated without the prior consent of Ocean Optics, Inc. in any form of binding or cover other than that in which it is published.

Trademarks

All products and services herein are the trademarks, service marks, registered trademarks or registered service marks of their respective owners.

Limit of Liability

Every effort has been made to make this manual as complete and as accurate as possible, but no warranty or fitness is implied. The information provided is on an "as is" basis. Ocean Optics shall have neither liability nor responsibility to any person or entity with respect to any loss or damages arising from the information contained in this document.

Ocean Optics. For more information visit www.oceanoptics.com.

Table of Contents

About This Manual	iv	Flame-NIR LEDs	10
Document Purpose and Intended Audience	iv	Change the Slit	10
Document Summary	iv	Accessories.....	11
Product-Related Documentation.....	v	Cables and Connectors	11
Document Version.....	v	DB15 Connector Cable (FLAME-CBL-DD4P-DB15P).....	12
Patent Pending Notice.....	v	PAK50 Connector Cable (FLAME-CBL-DD4P-PAK50P)	13
Warranty	v	Breakout Box (HR4-BREAKOUT)	13
ISO Certification	vi	Interchangeable Slits	13
Compliance	vi	Light Sources, Cuvette Holders and Other Accessories.....	14
 		Measurement Techniques – Typical Set-ups	14
Chapter 1: Introduction	1	Absorbance	14
Product Introduction	1	Common NIR Applications	15
Product Features.....	2	Reflectance & Transmission.....	15
Typical Applications.....	3	 	
Product Versions.....	5	Common NIR Reflectance and Transmission Applications	15
 		Irradiance.....	16
Chapter 2: Installation and Setup.....	7	Common Irradiance Applications	17
What's In the Box.....	7	 	
Flame-NIR Installation	7	Chapter 3: Flame-NIR Operation with OceanView	19
Software Installation.....	8	Overview	19
About OceanView.....	9	Launch OceanView	19
Hardware Setup.....	9	OceanView Main Screen.....	20
Hardware Feature.....	10	Connect the Flame-NIR in OceanView.....	21
		Set Acquisition Parameters	22
		Quick View and Device Response.....	23

Continuous and Single Acquisitions.....	23	Pixel Definition.....	46
Save Data	23	CCD Detector Reset Operation	46
Saved Data Panel.....	24	Timing Signals	47
Projects and Methods.....	25	Strobe Signals.....	47
Spectroscopy Application Wizards	25	Single Strobe.....	47
Dark and Reference Measurements	25	Continuous Strobe.....	47
Schematic View	26	Synchronizing Strobe Events	47
Chapter 4: Troubleshooting	29	External Triggering.....	48
Overview	29	Triggering Modes.....	48
Microsoft Windows Operating Systems.....	30	Normal	48
Apple Mac OSX Operating Systems.....	31	External Synchronous Trigger Mode.....	49
Linux Operating Systems.....	31	External Hardware Level Trigger Mode.....	50
Chapter 5: How the Flame-NIR Spectrometer Works	33	External Hardware Edge Trigger Mode.....	51
Overview	33	Chapter 7: Calibration.....	53
Chapter 6: Technical Specifications	37	Overview	53
Mechanical Diagram	40	Wavelength Calibration	53
Electrical Pinout.....	41	About Wavelength Calibration	53
DD4 Accessory Connector Pinout Diagram	41	Calibrating the Spectrometer Wavelength.....	54
SPI.....	44	Preparing for Calibration.....	54
I2C.....	44	Calibrating the Wavelength of the Spectrometer.....	54
Performance Charts for Flame-NIR	44	Irradiance Calibrations	57
Detector Overview	45	Chapter 8: Firmware and Advanced Communications	59
Detector Description	45	Flame-NIR Firmware.....	59
Detector Charge Capacity.....	45	Hardware Description	59
ii.....		USB Information	59
2 2 5 - 2 0 0 0 0 - 0 0 0 - 1 1 - 2 0 1 6 0 4		Instruction Set.....	59
Signal Averaging.....	46	Command Syntax	59
Internal Operation.....	46	USB Command Summary	60

USB Command Descriptions 61

Flame-NIR Serial Port Interface Communications and Control Information.... 73

Hardware Description 73

Instruction Set..... 73

Command Descriptions 75

Index 85

Copyright © 2026 Ocean Optics.
All rights reserved. No part of this publication may be reproduced, stored in a retrieval system, or transmitted, by any means, electronic, mechanical, photocopying, recording, or otherwise, without written permission from Ocean Optics.

This manual is sold as part of an order and subject to the condition that it shall not, by way of trade or otherwise, be lent, re-sold, hired out or otherwise circulated without the prior consent of Ocean Optics, Inc. in any form of binding or cover other than that in which it is published.

Trademarks
All products and services herein are the trademarks, service marks, registered trademarks or registered service marks of their respective owners.

Limit of Liability
Every effort has been made to make this manual as complete and as accurate as possible, but no warranty or fitness is implied. The information provided is on an "as is" basis. Ocean Optics shall have neither liability nor responsibility to any person or entity with respect to any loss or damages arising from the information contained in this document.

About This Manual

Document Purpose and Intended Audience

Thank you for choosing Ocean Optics! We hope that you'll be delighted with your decision. This document provides the users of Flame-NIR+ Spectrometers with instructions for setting up, calibrating and performing experiments with their spectrometer. It also contains detailed technical specifications and information about firmware and hardware integration. If you can't find what you're looking for in this document, please do not hesitate to contact us at techsupport@oceanoptics.com or via www.oceanoptics.com.

Document Summary

Chapter	Description
Chapter 1: Introduction	Introduces the product features. Contains descriptive information about the Flame-NIR Spectrometer. It also provides a list of system requirements, typical applications, and product versions.
Chapter 2: Installation and Setup	Provides installation instructions, including how to set up the Flame-NIR with OceanView. Also includes package contents and typical set-ups for different measurement techniques.
Chapter 3: Flame-NIR Operation with OceanView	Describes how to use the Flame-NIR with OceanView software, including how to connect, acquire, save and other basic features.
Chapter 4: Troubleshooting	Contains recommended steps to isolate and correct common problems.
Chapter 5: How the Flame-NIR Spectrometer Works	Describes how the Flame-NIR operates, illustrating the various parts and functions.
Chapter 6: Technical Specifications	Contains technical specifications and connector pinouts for the Flame-NIR Spectrometer.
Chapter 7: Calibration	Provides instructions for calibrating the Flame-NIR Spectrometer.
Chapter 8: Firmware and Advanced Communications	Contains a description of firmware commands.

Patent Pending Notice

The Flame-NIR spectrometer is covered by Patents Pending. Any violation of Ocean Optics intellectual property will be prosecuted.

Warranty

Our 3-Year Warranty covers Ocean Optics miniature fiber-optics spectrometers, spectral sensors, light sources and sampling accessories – regardless of the application – from manufacturing defects. It also covers fibers and probes for a full 12 months: <http://oceanoptics.com/services/exclusive-3-year-warranty/>.

This comprehensive warranty ensures you of the highest level of craftsmanship and reliability for years to come. No other manufacturer offers

such a solid guarantee of quality and reliability. The Ocean Optics 3-Year Warranty applies to Ocean Optics equipment (excluding OEM configurations) purchased on or after July 1, 2010. The warranty covers parts and labor needed to repair manufacturing defects that occur during the warranty period. We also will cover the costs of shipping warranty-related repairs from our customers to Ocean Optics and from us to our customers.

ISO Certification

Ocean Optics, the industry leader in miniature photonics, has been certified for ISO 9001:2008 certification applicable to the design and manufacture of electro-optical equipment since 2009.



Warning

This is a Class A product. In a domestic environment this product may cause radio interference in which case the user may be required to take adequate measures.



FCC Compliance

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which the user will be required to correct the interference at his own expense.



Warning

The authority to operate this equipment is conditioned by the requirement that no modifications will be made to the equipment unless the changes or modifications are expressly approved by the manufacturer.

Chapter 1: Introduction

Product Introduction

Combining the small size of the Flame optical bench with a new uncooled InGaAs detector, the Flame-NIR opens a new frontier in NIR spectroscopy. Without thermoelectric cooling, the Flame-NIR's power consumption is very low and makes the Flame-NIR ideal for integration into handheld and portable systems.

The Flame-NIR benefits from all the advantages of the Flame product line, including interchangeable slits and low unit-to-unit variability. With a preconfigured range from 970-1700nm, Flame-NIR is an ideal choice for many NIR applications including moisture measurement, grain and feed quality, measurement of fats and oils and pharmaceutical ingredients blending.



Flame-NIR+ Spectrometer

NIR spectroscopy is a powerful method for identification and classification as well as quantification and characterization of various parameters across a vast range of samples. NIR spectroscopy is fast, non-destructive and typically requires zero sample preparation. When combined with Ocean Optics extensive range of light sources and sampling accessories for reflectance, irradiance or transmission the Flame-NIR+ is ideal for many NIR applications (see Typical Applications).

Product Features

- Compact yet rugged. The small size of the Flame-NIR+, just 89.3 x 63.3 x 31.9 mm, makes it ideal for integrating into portable and handheld systems. With no moving parts the Flame is ideal for use in demanding environments.
- Sensitivity. Despite the compact footprint, the Crossed Czerny Turner design and InGaAs detector deliver superior sensitivity compared to other miniature NIR designs. Measure more samples faster using the Flame-NIR+.
- Ease of use. Just plug in the micro-USB connector and start measuring. Design experiments and projects in OceanView software. Modular accessories make it easy to change between reflectance, transmission/absorbance or irradiance while optimizing for specific samples and applications.
- User interchangeable slits. Take control with interchangeable slits that allow the user to vary the throughput and resolution in minutes. Swap slits between SMA and FC type connectors to work with your application.

- Low unit to unit variation. Cutting edge manufacturing techniques have reduced unit to unit variance in almost every aspect compared to previous generations of the spectrometer, including uncalibrated sensitivity.
- Low power consumption. Drawing <250mA at 5V, the Flame-NIR+ can be powered almost anywhere including in OEM portable and handheld devices and industrial process control environments.
- Indicator LEDs show the power and data transfer status of the spectrometer at all times.
- Choice of USB or RS-232 serial communications supported by drivers and software that make it easy to integrate into almost any system.
- 8 GPIO pins and the ability to provide power to external devices through the 40-pin connector.
- Purchase with complete peace of mind with Ocean Optics industry leading 3-year guarantee.

KeyFeature	BestFor...	ExampleApplications
User-Interchangeable Slits	Users who wish to vary resolution and throughput during measurements, or switch from absorbance to fluorescence in minute	Life science and other labs using a wide variety of biological samples
Indicator LEDs	Convenient visual reference for spectrometer operation and status	Teaching and general lab use
Thermal Stability	Applications that require repeatable results in industrial and other environments with varying temperatures and conditions	LED binning and light metrology, process monitoring
Reduced Unit-to-Unit Variation	OEM applications such as uncalibrated sensitivity and other measurement needs where users benefit from low variance	OEM manufacturers of analytical instrumentation
Configurability	Optimizing your spectrometer for application-specific requirements; adjusting range, throughput and resolution and adding features such as filters as required	Laser characterization, low signal fluorescence and optimizing for specific absorbance bands
Plug & Play Operation	Users that want the convenience of simple, fast operation via the micro USB connection; lets users take the measurement to the sample	Remote sensing measurements in the field, including air and water quality monitoring and solar irradiance
Ease of Integration	OEMs and developers who need to integrate a spectrometer as part of a system via USB or RS-232	Engineering labs, developers, OEM manufacturers; works with LabVIEW and other design platforms

Installation and Setup

Typical Applications

<i>Application Area</i>	<i>Examples</i>
Food & Agriculture	• Freshness and sweetness of fruits and vegetables • Fat and protein content of meat • Grain quality parameters including moisture, starch and from farm to table • Identification and classification of seeds, grains and pulses
Life Sciences & Pharmaceuticals	• Identification of Pharmaceutical Ingredients • Portable diagnostic equipment for blood glucose and blood flow • API blending in Pharmaceutical manufacturing
Light & Laser	• Measure Laser center wavelength and stability • Monitor NIR irradiance and background in and out of the lab
Chemical & Commodity Manufacturing	• Create custom chemometric models to measure quality parameters in hydrocarbons extraction and processing • Identification of plastics of various type and color in recycling

Specifications Summary

Specification	Flame-NIR+
SPECTROSCOPIC	
Optical resolution	11 μ m (25 μ m slit)
Signal-to-noise ratio	6000:1
Integration time	1 ms – 3 seconds*
Corrected linearity	99.8%
Range	970 nm - 1700 nm
ELECTRONIC	
A/D resolution	16 bit
Power requirement (spectrometer functions)	250 mA @ +5 VDC
Inputs/Outputs	8 x digital user programmable GPIOs
Trigger modes	4 modes (see Triggering Modes)
Strobe functions	Yes
Gated delay feature	Yes
Connectors	Micro-USB and JAEDD440-pin connector
DETECTOR	
Detector	Hamamatsu G8160-03 InGaAs linear image sensor
Detector range	900 – 1700 nm
Pixels	128 pixels
Invalid Pixels	Minimum: 0 Maximum: 2
PHYSICAL	
Dimensions	88.9 mm x 63.5 mm x 31.9 mm (excluding feet)
Weight	265 g
* Because the baseline will rise rapidly with temperature and extended integration times Ocean Optics recommends integration time be kept below 3 seconds when using the Flame-NIR.	

Product Versions

Ocean Optics offers both preconfigured units as well as custom-configured units, enabling you to order a customized spectrometer optimized for your application. You can tell which kind of spectrometer you have by looking at the product code, located on the bottom of your spectrometer.

You can find more information about the various components and possible configuration options in Chapter 5: How the Flame-NIR+ Spectrometer Works.

Product Code Taxonomy

Spectrometer Family?	Detector Type?	Preconfigured or Custom?
FLAME-	NIR (Hamamatsu G8160-03)	-INTSMA25 (preconfigured)
		-XXX(custom)

Preconfigured Models

Flame-NIR+Preconfigured Models (Hamamatsu G8160-03 Detector)								
	Model	Range (nm)	Std Slit (μm)	Resolution std slit (nm)	Grating #	Mirror Type	Filters & Detector Options*	Lens
General Purpose	Flame-NIR-INTSMA25	970-1700nm	25	11nm	NIR33	Au	OF1-830g	no

For more information and specifications on preconfigured models, see www.oceanoptics.com.

Chapter 2: Installation and Setup

What's In the Box

- **Packing List**

The packing list is inside a plastic bag attached to the outside of the shipment box (the invoice arrives separately). It lists all items in the order, including customized components in the spectrometer (such as the grating, detector collection lens, and slit). The packing list also includes the shipping and billing addresses, as well as any items on back order.

- **Flame-NIR+ Spectrometer**

Your Flame-NIR+ spectrometer arrives pre-calibrated and ready to plug and play.

- **Micro USB Cable**

Use this cable (CBL-MICROTOA-USB) to connect your spectrometer to a computer running on a Windows, Mac or Linux operating system.

- **Wavelength Calibration Data Sheet**

Each spectrometer is shipped with a Wavelength Calibration Data Sheet that contains information unique to your spectrometer. OceanView reads this calibration data from your spectrometer when it interfaces to a computer via the USB port.

Note: Please save the Wavelength Calibration Data Sheet for future reference.

Flame-NIR+ Installation

The following procedure provides general instructions for getting your new Flame spectrometer up and running.

CAUTION




Be sure to install the software BEFORE connecting the spectrometer to your PC. The software installs the drivers required for spectrometer installation. If you do not install the software first, the system will not properly recognize the spectrometer.

If you have already connected the Flame to a computer running on a Windows platform prior to installing the operating software, consult Chapter 4: Troubleshooting for information on correcting a corrupt Flame installation.

Procedure

Use the following procedure to get your system up and running quickly. See Hardware Setup for more detailed information.

1. Install your spectroscopy operating software by following the installation wizard prompts. See Software Installation for more in-depth information.
2. Locate the micro USB cable provided with the spectrometer. Connect the spectrometer to the USB port on your computer using this cable.

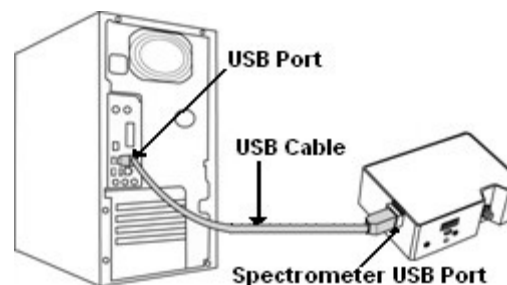

3. Take measurements. Refer to your spectrometer operating software manual for information on the software user interface.

Software Installation

Use OceanView version 1.5.2 and above for Flame-NIR+. Flame-NIR+ cannot be used with SpectraSuite. You can use OceanView on the following operating systems.

CAUTION



Be sure that you download the correct software package for your computer version (32 or 64-bit). See the Frequently Asked Questions in Chapter 4: Troubleshooting for more information on determining your computer version.

Software OS	Windows							Apple	Linux
	2000	XP	Vista	7	8	8.1	10		
OceanView		✓	✓	✓	✓	✓	✓	OSX Version 10.5 or later on Intel processor	Any version released for x86 or amd-64 platform since 2010


About OceanView

OceanView is the latest generation of operating software for all Ocean Optics spectral devices. It is a Java-based spectroscopy software platform that operates on Windows, Macintosh and Linux operating systems. The software can control any Ocean Optics USB device.

OceanView is a user-customizable, advanced acquisition and display program that provides a real-time interface to a variety of signal-processing functions. With OceanView, you have the ability to perform spectroscopic measurements (such as absorbance, reflectance, and emission), control all system parameters, collect and display data in real time, and perform reference monitoring and time acquisition experiments. Consult the OceanView manual for hardware requirements when using OceanView (see Product-Related Documentation).

Hardware Setup

The Flame-NIR Spectrometer connects to a computer via the USB port or serial port. When connected through a USB 2.0 or 1.1 port, the spectrometer draws power from the host computer, eliminating the need for an external power supply. The Flame-NIR, like all Ocean Optics USB devices, can be controlled by our OceanView software (see About OceanView).

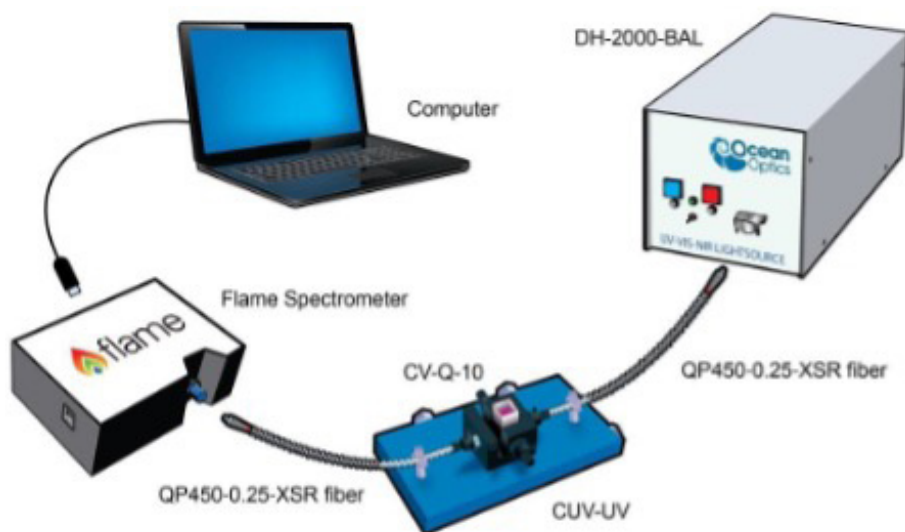

Ocean Optics Flame-NIR+ Fiber Optic Spectrometer Typical Set-up

Follow the steps below to connect the Flame-NIR+ to a computer via the USB port:

1. Install the spectrometer operating software on the destination computer.
2. Locate the USB cable (CBL-MICROTOA-USB) provided with the Flame-NIR+.
3. Insert the micro connector end of the cable into the side of the Flame-NIR+ and the larger end of the cable into the USB port of the computer.
4. Connect any spectroscopy accessories. To find operating instructions for Flame-NIR+-compatible products (such as light sources, sampling chambers, and probes), consult the Ocean Optics website.
5. Attach the fiber to the fiber optic connector on the spectrometer.

If you installed the spectrometer operating software prior to connecting the Flame-NIR, the software automatically installs the Flame-NIR drivers. If the drivers do not successfully install (or if you connected the Flame-NIR to the computer before installing the software), consult Chapter 4: Troubleshooting.

Hardware Features

Flame-NIR LEDs

The Flame-NIR features two indicator lights that operate as shown below:

LED	Steady	Flashing
red	Unit is on	N/A
green	Unit is ready	Unit is acquiring data

Note that LEDs can be turned off in OceanView or by using a firmware command.

Change the Slit

The Flame-NIR allows you to change your spectrometer's slit to match your application requirements. Note that a filter must be ordered for each slit (if your application requires a filter). There's no need to calibrate your spectrometer when changing the slit, just install and start measuring!

• Procedure

1. Find the SMA connector. If a fiber is attached, remove it.



1. Use the Allen key to remove the 2 the screws attaching the slit to the spectrometer.
2. Pull the slit out of the spectrometer.
3. Put the new INTSMA slit connector into the spectrometer; with the key of the connector on the left side.
- 4.. Put the new INTSMA slit connector into the spectrometer; with the key of the connector on the left side.
5. Install the 2 screws again. Use the Allen key to tighten the screws carefully (do not over-tighten).
6. If necessary, connect the fiber again.




Cables and Connectors

Cables are available to connect your Flame-NIR+ Spectrometer to accessories:

- Accessory cable for light sources and other accessories (DB15 Connector Cable)
- Accessory cable for HR-type connector (PAK50 Connector Cable)
- Breakout board (DD4 -BREAKOUT-BOARD) This breaks out the 40 pin DD4 connector to a set of header pins that can be used to wire each pin, as required.

DB15 Connector Cable (*FLAME-CBL-DD4P-DB15P*)

This cable connects the Flame-NIR+ to existing Ocean Optics accessories that use a DB-15HD connector. These include the PX-2, LLS and HL-2000-FHSA light sources.

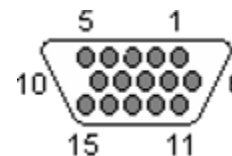

DD4-DB15PinConnections

Accessories

Ocean Optics provides a range of standard cables and accessories that connect the Flame-NIR+ to our large range of sampling and light source accessories. Items specifically designed for the Flame-NIR+ are described here; they are not provided with the Flame-NIR+ spectrometer and must be purchased separately.

15PIN	DD4	Name
1	4	SingleStrobe
2	3	ContinuousStrobe
3	40	Vusb
4	2	ExternalTriggerIn
5	2	ExternalTriggerIn
6	7	GPIO1
7	NC	Reserved(NC)
8	2	ExternalTriggerIn
9	8	GPIO2
10	1	Ground
11	22	I2CSDA
12	21	I2CSCL
13	5	LampEnable
14	NC	Reserved(NC)
15	11	GPIO4

PPAK50 Connector Cable (FLAME-CBL-DD4P-PAK50P)

This cable connects the Flame-NIR+ to the Breakout Box.

DD4P to PAK50P 30-Pin Connections

PAK50P	DD4	Name	PAK50P	DD4	Name
1	25	RS232 Rx	16	11	GPIO 4
2	24	RS232 Tx	17	4	Single Strobe
3	8	GPIO 2	18	12	GPIO 5
4	40	VUSB	19	16	SPI CLK
5	1	GND	20	3	Continuous Strobe
6	21	I2C SCL	21	18	SPI CS
7	6	GPIO 0	22	13	GPIO 6
8	22	I2C SDA	23	28	Reserved (do not connect)
9	7	GPIO 1	24	NC	Reserved (NC)
10	2	External Trigger In	25	5	Lamp Enable
11	9	GPIO 3	26	14	GPIO 7
12	40	VUSB	27	10	GND
13	17	SPI MOSI	28	NC	Reserved (NC)
14	40	VUSB	29	15	GND
15	19	SPI MISO	30	NC	Reserved (NC)

For the DD4, connect shield to connector case, if shield is included.

Breakout Box (HR4-BREAKOUT)

The Breakout Box is a passive module that separates the signals from the Flame-NIR+'s DD4 40-pin connector to an array of standard connectors and headers, enabling functionality with a wide range of accessories. In addition to the accessory connector, the breakout box features a circuit board based on a neutral breadboard pattern that allows custom circuitry to be prototyped on the board itself. See Product-Related Documentation to access the manual for the Breakout Box.

Interchangeable Slits

The Flame-NIR+ offers the capability of changing the slit size to match your measurement needs. You can order additional replacement slits either individually or as a kit (in various widths from 5 μm to 200 μm). See Change the Slit and Chapter 5: How the Flame-NIR+ Spectrometer Works for more information.

Light Sources, Cuvette Holders and Other Accessories

Ocean Optics supplies a large range of accessories for use with our spectrometers.

- Fibers
- Light Sources
- Integrated Sampling Systems
- Cuvettes, including microfluidic cuvettes
- Filter Holders & Filters including Low Pass, Band Pass and High Pass


Measurement Techniques – Typical Set-ups

The Flame-NIR+, in conjunction with Ocean Optics light sources and sampling accessories, can be used for many different measurement techniques. One of the key advantages of modular fiber optic spectroscopy is that you can change components of the system without having to buy a whole new system. Here, we show a range of typical UV-Vis set ups for basic spectroscopy techniques. You'll find lots more information about measurement techniques at www.oceanoptics.com/measurementtechnique.

Absorbance



Absorbance is typically a relative measurement, comparing the spectrum from the sample to that of a reference. Absorbance is commonly used for concentration measurements and for identifying components in mixtures. The absorbance measurement scales the response logarithmically. Connect the Flame-NIR+ to our cuvette accessories via the SMA Adaptor accessory to take a liquid sample Absorbance measurement, or mount it directly against the sample with a light source on the opposite side for solid sampling.

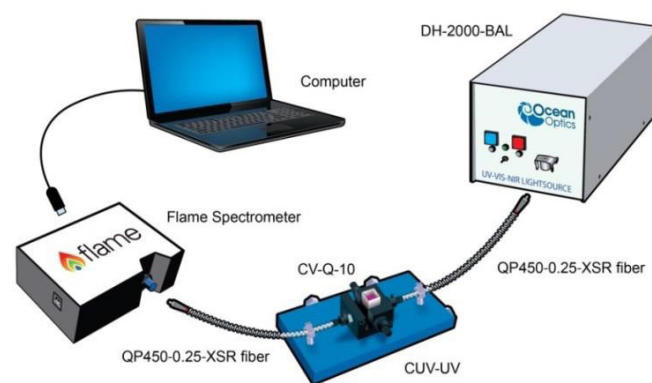

Typical Absorbance Set Up


Common NIR Applications

- Quantification of blood oxygen levels using a ratio of absorbances
- Concentration of solutions & gaseous samples such as CO₂
Absorbance scaling is used in many chemometric applications of NIR spectroscopy to highlight small shifts in spectral response

Reflectance & Transmission



Reflectance spectroscopy compares the relative level of light reflected off a sample compared with a reference (given as a percentage of the reference spectrum at each wavelength). A reflectance standard is used to set the reference level of 100%. Transmission is similar but compares the light transmitted through a sample, relative to a reference, rather than reflected off it. Typically Reflectance uses a fiber optic probe attached to a light source and a spectrometer, but measurements can be done easily with the Flame-NIR+, both in free-space or with the SMA Adaptor accessory. Transmission setups are usually the same as Absorbance setups.

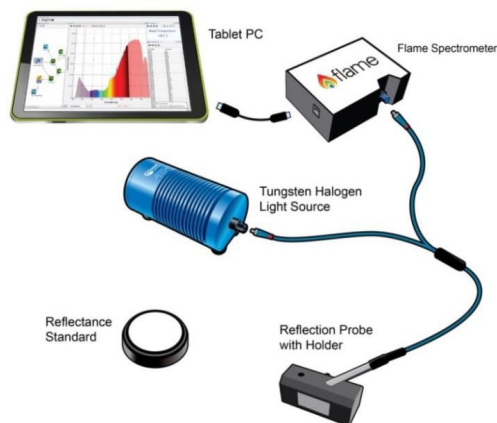

A Reflectance Set Up with Probe, Reflectance Standard and Probe Holder

Common NIR Reflectance and Transmission Applications


- Freshness and sweetness of fruits and vegetables
- Fat and protein content of meat
- Grain quality parameters including moisture, starch and from farm to table
- Identification and classification of seeds, grains and pulses
- Identification of pharmaceutical ingredients
- Portable diagnostic equipment for blood glucose and blood flow
- API blending in pharmaceutical manufacturing
- Thin film and semiconductor metrology

Irradiance



Irradiance is the technique of measuring the total energy of light at a given wavelength, either relative to the spectral output of a known source (relative irradiance) or in absolute units of power or energy (absolute irradiance). This is used widely in light metrology, color measurement and environmental science. Absolute irradiance measurements require an irradiance-calibrated spectral device. This can be done in the factory for some configurations or by using a calibration lamp in the lab or field. Every time a set-up is changed, the device used must be recalibrated.


For a more detailed explanation of how to calibrate a device see Chapter 6: Calibration, or contact Ocean Optics. Use the front-mounted diffuser and SMA adapter accessories for irradiance measurements with the Flame-NIR+.

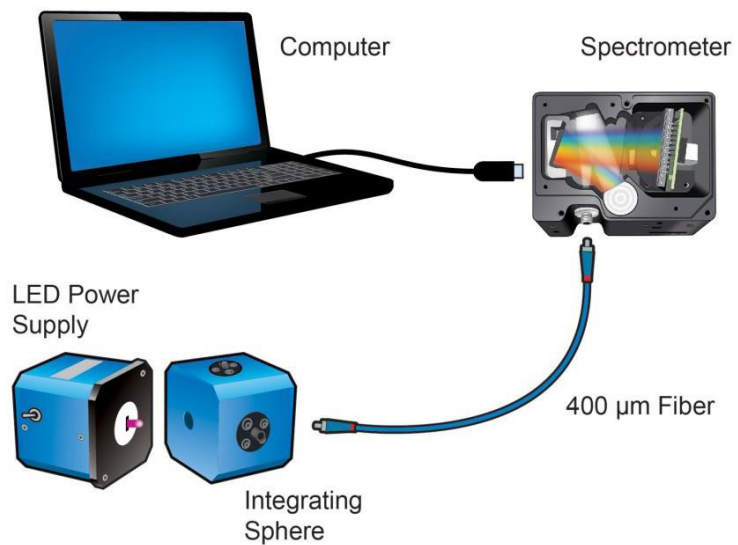

Typical Relative Irradiance Set-up for Measuring Light Power Output of an LED Using and Integrating Sphere

Common Irradiance Applications

- Measuring the radiant output of lamps and LEDs
- Measuring laser center wavelength and stability
- Monitoring NIR irradiance and background in and out of the lab

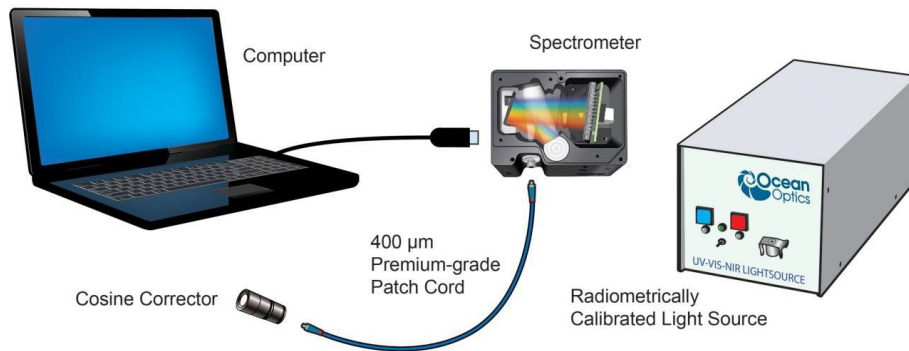

Typical Set-up for an Absolute Irradiance Measurement Using Field Calibration with a Calibrated Light Source

Chapter 3: Flame-NIR+ Operation with OceanView

Overview

The following information enables you to perform the basics of acquiring and saving data with your Flame-NIR+ Spectrometer and OceanView software. More detailed information about OceanView is in the OceanView Manual.

Launch OceanView

Once you have installed your software and connected your spectrometer, you are ready to display your measurement data using OceanView. Launching OceanView differs, depending on your operating system and where you have placed your OceanView program files.

For PCs running Microsoft Windows, the default location is Start | Programs | Ocean Optics | OceanView | OceanView.

For Mac computers, the default location is the Applications folder.

When you first start OceanView, the Welcome Screen appears.



The OceanView Welcome Screen (Version 1.5)


Quick View - Displays the spectrum in Quick View mode showing raw, unprocessed data. This is uncorrected for instrument response vs. wavelength. Quick View shows you a live shot of what the Flame-NIR+ is "seeing." From Quick View you can launch application wizards or construct your own method.

Load a Saved Project - Loads a previously saved project. Click Restore Last Session to reload the schematic and views as they were when the software was last closed.

Spectroscopy Application Wizards – Use this function to set up a measurement using simple step-by-step wizards. A large range of applications is available.

OceanView Main Screen

No matter what route you take on start up, you will soon end up on the OceanView main screen. This is where you can set and view acquisitions, save and load data and save projects.

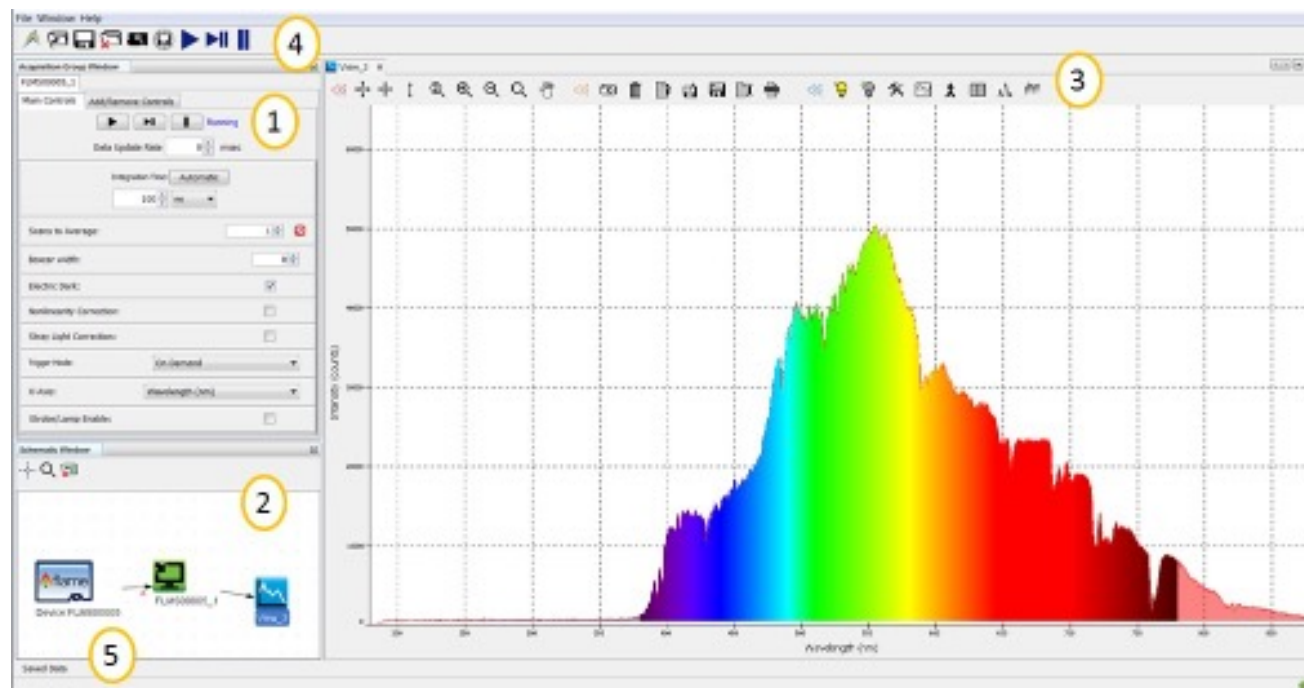

1. Acquisition Group Window	Use to set acquisition parameters such as integration time. Controls the spectrometer acquisition.
2. Schematic View	Schematic view graphically displays the flow of information from the spectrometer to the view. Use nodes to mathematically modify the data to create processed measurements (methods). This function is extremely flexible and incredibly powerful.
3. View Display	Display your data, view, save and display controls, as well as other features such as peak finder and quick dark & reference.
4. Global Controls	Control all spectrometers synchronously, save projects, and start a new application wizard.
5. Saved Data	Displays data saved in the active save file path. Preview data, store notes and load overlays directly to the active view. Click to open.


Connect the Flame-NIR+ in OceanView

The Flame-NIR+ should automatically appear when you start OceanView and should be acquiring with the default acquisition parameters. If you do not see a signal or the Flame-NIR+ icon on the schematic you may need to rescan for spectrometers.

Procedure

To rescan for attached devices,

1. Click on the Device Manager icon ().
2. Click Rescan. The spectrometer should automatically connect.

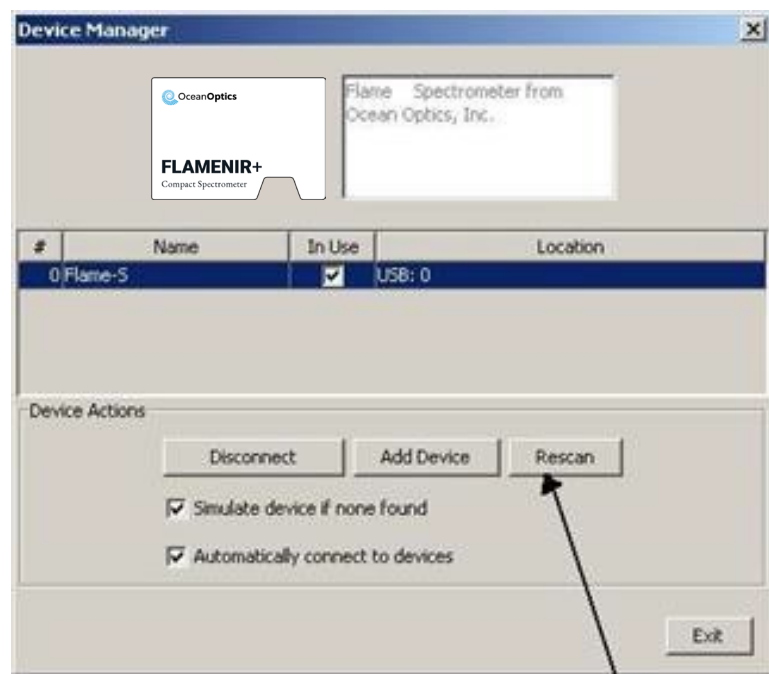

Rescan Button

Set Acquisition Parameters

Set Acquisition parameters in the Acquisition Group Window to control the spectrometer. This window may be minimized when you first start OceanView. You can either expand or open a new window from the menu (Window | Acquisition Group). An active acquisition is required for the Acquisition window to appear. Functions available to control in the Acquisition window include the following:

- **Integration Time** – Sets the integration time, the time over which the detector captures incident light. At the end of the integration time the accumulated signal is read from the detector by the electronics.
- **Averaging** – Signal, especially at low levels, is often significantly impacted by noise. Averaging several spectra together reduces the impact of noise and provides a cleaner result. However, at long integration times, averaging can increase the total time of a measurement significantly.
- **Boxcar** – Boxcar is a form of averaging across pixels. It applies a rolling average to multiple adjacent pixels to help smooth the spectral response and reduce the impact of noise.
- **Non Linearity Correction (on/off)** – Detectors do not have a completely linear response. As they approach saturation, typically their efficiency reduces.
- **Trigger Modes** – Sets triggering mode. For more information on triggering see External Triggering.
- **Strobe/ Lamp (on/off)** – Use this function to turn an attached light source on or off.
- **GPIO Controls** – Can be used to control compatible accessories or custom hardware. Can be set to three states, on, off and alternate.

Controls that appear in this window depend on the spectrometer model. You can add and remove acquisition controls from this window.

Quick View and Device Response

In **Quick View mode** (formerly **Scope mode**) the spectra that are displayed have an arbitrary y-axis given in counts. This is the raw signal from the detector and is proportional to the voltage induced by the light falling on the detector.

It is very important to realize that this is uncalibrated data and that a counts signal does not represent a particular power or energy from one wavelength to the next. Because the response of the detector is linear, twice the counts at a particular wavelength do indicate that the amount of light at that wavelength has “doubled” (relative to another wavelength). However, a small peak relative to a big peak does not indicate that there is less or more light at a particular wavelength relative to another in absolute terms. To understand the true relationship you need to do a relative measurement (including relative irradiance), or if you want a quantified result, an absolute irradiance measurement.

The relative efficiency of converting light to detector signal varies significantly across the range of a spectrometer. Many things impact this including the responsiveness of the detector (quantum efficiency) and the efficiency of optical components. Each configuration of a spectral device has a unique response curve, referred to as the instrument response.

Continuous and Single Acquisitions

There are two sets of controls for taking or pausing acquisitions. The set on the Acquisition group window allows you to control each device individually. The set on the top bar is a global control that will allow you to start and pause all devices currently attached.

	Acquire data continuously
	Take a single acquisition and then pause
	Pause all acquisitions.


Save Data

	Configure Saving, set saving parameters and file type, file directory and file naming convention. Once selected, the file directory will persist until changed.
	Start saving data. Turns red when save is active. If saving data continuously, click when red to stop saving. Will only activate saving for acquisitions attached to that particular view.
	Global Save. Activates all configured saves across all views. Use to save data from multiple devices at the same time.

By default OceanView will save data as a single “snapshot” acquisition. By configuring the save you can set the save behavior to fit your measurement needs, from single snap shots to a continuous stream of data over time.

Saved Data Panel

The saved data panel lets you see your data as it is saved and preview data. It also makes it simple to add overlays of saved data to your screen.

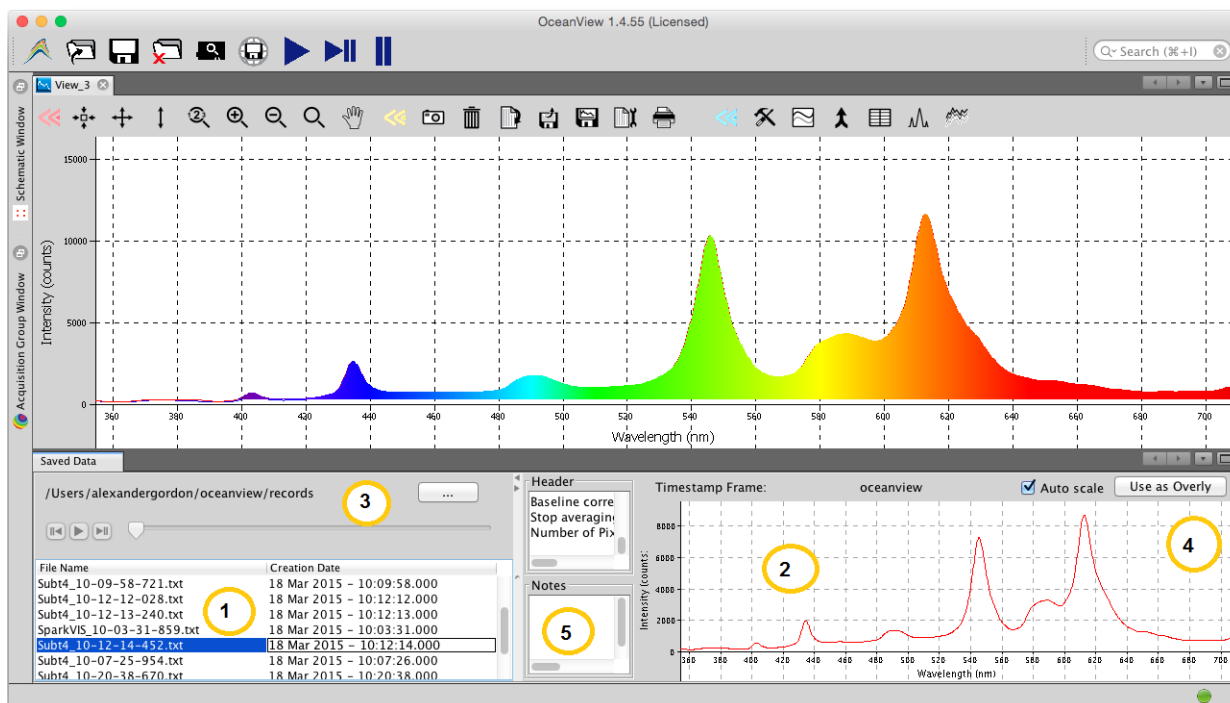

Saved Data Panel

1. Saved Files	List of saved files currently in the saved directory. Arrange by name or date.
2. Preview	Shows a preview of the saved spectra, time series or appended series saved data can be stepped through acquisition by acquisition using the controls above the saved files list.
3. File Path	Set the file directory.
4. Overlay	Set the previewed spectra as an overlay on the active view.
5. Notes	Enter notes about the saved spectra. Notes are saved with the same file name as a separate .tsv file. These can be viewed or edited with any text viewer such as notepad.


Projects and Methods

OceanView makes it easy to save and load projects and methods. We define a project as a measurement set up made with a particular spectral device. If the software cannot find the device, it will load this as a method and prompt the user to select a substitute device from those selected.

	Click to save a project. Alternatively select <i>File Save Project</i> from the menu. Saves all view and schematic parameters to a single ASCII file.
	Load a project or method.

Spectroscopy Application Wizards

	Click this button to set up a measurement using simple step by step wizards. A large range of applications is available.
---	--

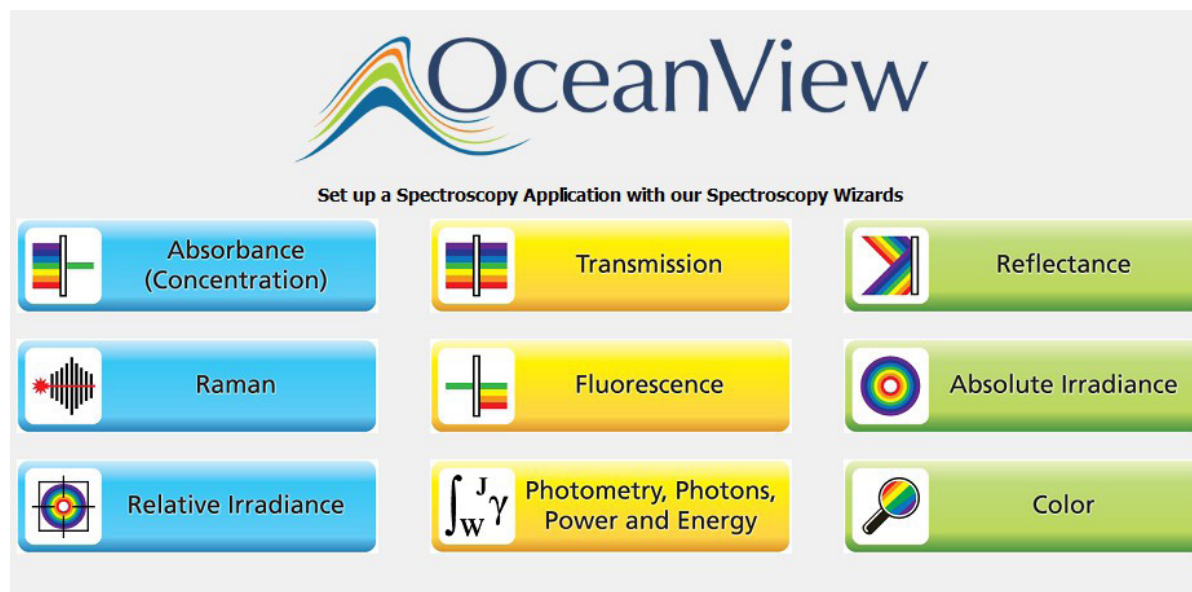

Application Wizard Window

Dark and Reference Measurements


Dark and reference measurements are commonly used in spectroscopy.

- **Dark Measurements** – subtract a background signal from the spectrum. This can be considered the removal of a constant error. Typically this is done when the light source is off to remove any background from the ambient environment, hence the name dark.
- **Reference Measurements** – make the signal relative to the reference. Consider this a normalization of the signal against a reference. Typically this is taken with a reference sample and the light source turned on. This lets you look at the relative spectral change compared to a reference sample.

Most often you will set up your measurement with the reference and dark through the application wizards. The wizards will prompt you to take your reference and darks. Alternatively you can use the quick dark and quick reference features. Once a dark and/or reference measurement has been set, you can update it with the controls on the top bar of the view.

	Quick Reference – click to take a reference and set up a new view. After clicking it will prompt the user to take a dark.
	Quick Dark – click to take a dark measurement and sets up a new Quick View minus dark view.
	Reference – click to update the stored reference measurement.
	Dark – click to update the stored dark measurement.

Schematic View

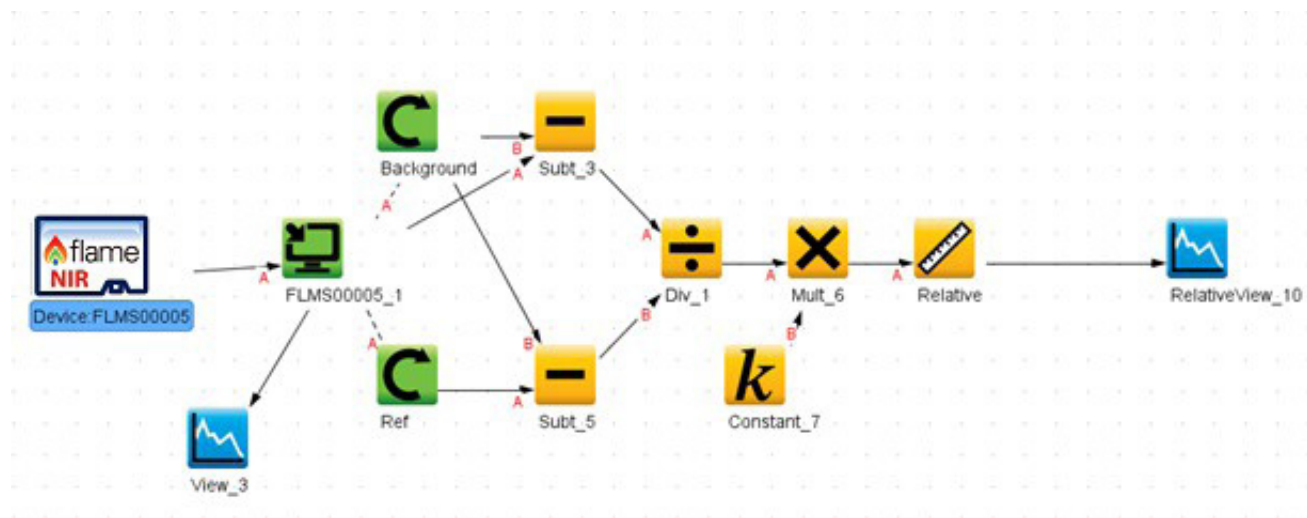

The schematic view is a graphical interface that allows you to move from device through to processed data. There are a few basic components to consider.

	Devices – Each spectrometer will appear as a separate device. Right click to open a menu that can generate an acquisition, control a TEC (if applicable) and add other device controls.
	Acquisitions – A spectrometer can output one acquisition per detector channel. Right click to open menu.
 Mult_6	Nodes – These are the building blocks of the schematic view. They are all various functions that take data in and provide an output. To make a node, right click on the schematic background. Each node can be configured by double clicking on the node. To join nodes press ctrl, click and drag (windows).
	Views – Are a type of window that displays data. To generate a new view right click on the schematic background.

More information about schematic view including detailed descriptions of the available nodes can be found in the *OceanView Installation and Operation Manual* and in the help section of the OceanView software.


Chapter 4: Troubleshooting

Overview

Sometimes things don't quite go to plan; hopefully you'll find some answers below. If not, don't hesitate to contact us and our Tech Support team will leap into action. Some typical questions are answered here. For more information, consult the FAQs on the Ocean Optics website at <http://oceanoptics.com/faq/>.

I connected the USB cable and started OceanView but I don't see my spectrometer attached.

Use the Rescan button in the Device Manager to rescan for attached devices.

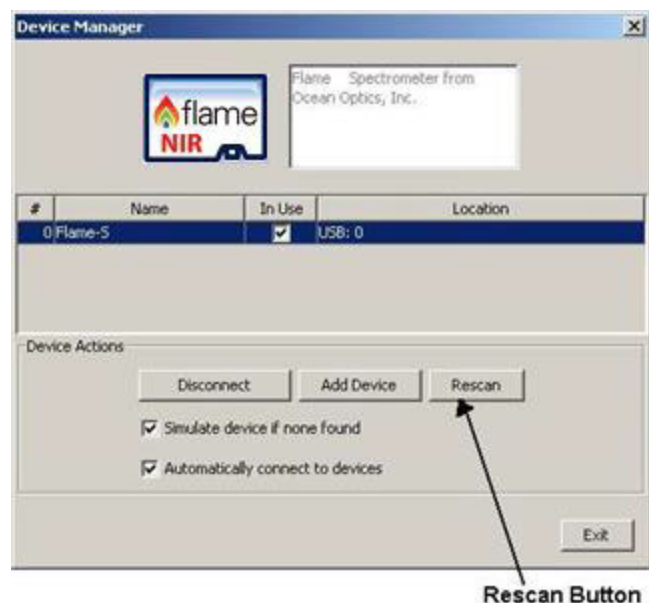

I am having trouble installing the drivers, what should I do?

Hardware device driver installation is usually seamless on Microsoft Windows operating systems and should happen in the background when you connect your spectrometer to a computer with the software installed. However, some Windows systems require a bit more care when connecting your spectrometer for the first time.

If your spectrometer is not recognized by OceanView on your computer, you need to manually install the spectrometer drivers. See your OceanView manual for this procedure. Also consult the Correcting Device Driver Issues document (see Product-Related Documentation).

How do I know my spectrometer has power?

The red LED on the spectrometer should be on steadily if the unit is receiving power.

How do I know my spectrometer is transmitting data?

The green LED on the spectrometer flashes when transmitting data.

I'm installing OceanView but I need a product key; where can I find this?

The product key was sent to the contact e-mail on the sales order when you purchased your OceanView license. Contact info@oceanoptics.com for more information. You'll need your sales order number, quotation number, the serial number of the spectrometer that was purchased with the software, and, if known, the e-mail address under which your product key was created to recover your key.

I connected the Flame-NIR+ to the computer before installing my spectroscopy operating software to install the drivers. What do I do now?

The steps to take to resolve this issue differ, depending on your computer's operating system.

Microsoft Windows Operating Systems

If you connected your Ocean Optics Flame-NIR+ device to the computer prior to installing your Ocean Optics software application on a Windows platform, you may encounter installation issues that you must correct before your Ocean Optics device will operate properly.

Follow the applicable steps below to remove the incorrectly installed device, device driver, and installation files.

Remove the Unknown Device from Windows Device Manager

- **Procedure**

1. Open Windows Device Manager. Consult the Windows operating instructions for your computer for directions, if needed.
2. Locate the **Universal Serial Bus Devices** option and expand the **Universal Serial Bus Devices** selection by clicking on the "+" sign to the immediate left.

Note: Improperly installed USB devices can also appear under the Universal Serial Bus Controller option. Be sure to check this location if you cannot locate the unknown device.

3. Locate the unknown device (marked with a large question mark). Right-click on the **Unknown Device** listing and select the **Uninstall** or **Remove** option.
4. Click the **OK** button to continue. A warning box appears confirming the removal of the Unknown Device. Click the OK button to confirm the device removal.
5. Disconnect the Flame-NIR+ from your computer.
6. Replug the spectrometer into your computer.

The system should now be able to locate and install the correct drivers for the USB device.

Apple Mac OSX Operating Systems

Since there are no device files for the Flame-NIR+ Spectrometer in a Mac operating system, you should not encounter any problems if you installed the spectrometer before the spectrometer operating software.

Linux Operating Systems

For Linux operating systems, all you need to do is install the spectrometer operating software, then unplug and replug in the spectrometer. Technically, the driver files for Linux simply give nonprivileged users permission to use newly connected hardware. There isn't any long-term harm to plugging in the device before installing the software.

How do I determine whether my Windows computer is 32-bit or 64-bit?

Errors can occur if you download the wrong version of software (for 32-bit or 64-bit computers). To verify your computer version for most Windows computers, go to the Properties window (under Computer or My Computer). If no version is listed, then your system is a 32-bit. For more information see the Microsoft Frequently Asked Questions at <http://windows.microsoft.com/en-us/windows/32-bit-and-64-bit-windows#1TC=windows-7>.

How do I check the configuration of my spectrometer?

Check the label on the bottom of your spectrometer. You can also check your configuration using your spectrometer operating software.

In OceanView: Open the Schematic window and double click the spectrometer icon.

Product Upgrades, Repairs and Servicing

Occasionally, you may find that you need Ocean Optics to make a change or an upgrade to your system. To facilitate these changes, you must first contact Customer Support and obtain a Return Merchandise Authorization (RMA) number. Please contact Ocean Optics for specific instructions when returning a product.

Repairs

Sometimes accidents happen! If you need to return your Ocean Optics Product for repair, here is what to do:

- **Procedure**
 1. Contact us to speak to an Ocean Optics representative about the problem. If it is determined that the product must be returned, the representative will issue an RMA number.
 2. Package your product, ideally in the original packaging, and return it to Ocean Optics, along with the RMA number that you received.

Note: For RMA returns under warranty we will organize and pay for shipping both ways. For accidental damage, you only pay to have the product delivered to your closest Ocean Optics or OOI Distributor Office.

Upon careful examination, we'll advise you with an estimate. When your product is ready, it will be returned to you.

Servicing


To keep your instrument in tip top shape we recommend yearly wavelength recalibration. You can do this yourself if you have appropriate tools or we can do this for you. Contact your local representative to find out more about service availability and cost. We offer the following services:

- Wavelength Calibration
- Absolute Irradiance Calibrations

Chapter 5: How the Flame-NIR+ Spectrometer Works

Overview

This section provides an overview of the Flame-NIR+ spectrometer and how it works from light entering the slit through to the transmission of the spectrum over USB. It also provides an overview of all the different possible configurations that are possible, designed to help you optimize your spectrometer for specific applications.

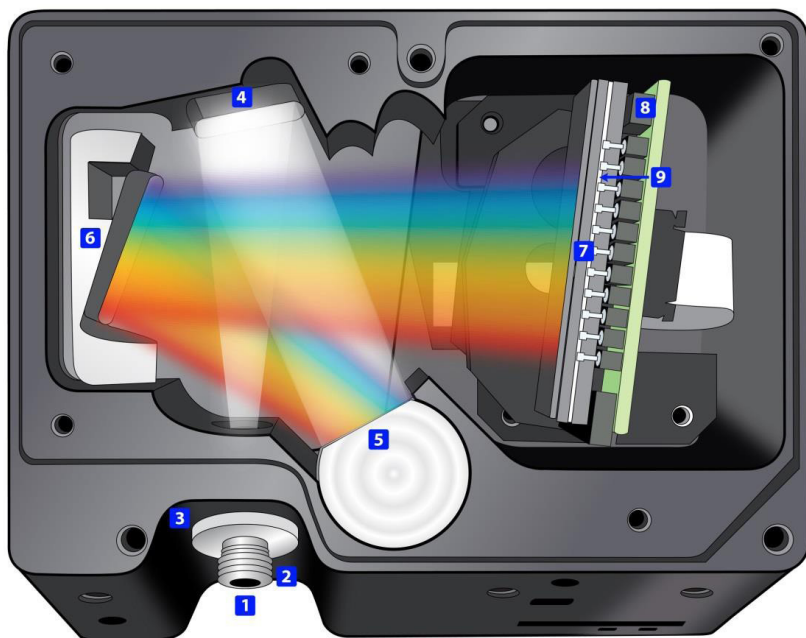

Flame-NIR+ Open Bench

1. **Fiber Optic Connector:** Light from a fiber enters the optical bench through the SMA 905 Connector. The SMA 905 bulkhead provides a precise location for the end of the optical fiber, slit, absorbing filter and fiber clad mode aperture. While we supply SMA connectors as standard, FC connectors are also available. See #2 for available options.
2. **Interchangeable Slit:** Light passes through the installed slit, which acts as the entrance aperture. Slits come in various widths from 5 μm to 200 μm . The slit is fixed in the SMA 905 bulkhead to sit against the end of a fiber. Smaller slit sizes achieve the best optical resolution while larger slits have higher light throughput. Slit size is labeled as shown.
3. **LongPass Absorbing Filter (optional):** If selected, an absorbing filter is installed between the slit and the aperture in the SMA 905 bulkhead. The filter is used to limit bandwidth of light entering spectrometer or to balance color. Filters are installed permanently. A filter is for a specific slit. If you anticipate needing the filter with multiple slit sizes, then you must specify this at the time you order. You will know which filter is installed in each slit because of the color-coded dots on the outside as shown in the figure and described in the table in the next page.



Item Code	Description	Dot 1	Dot 2
OF1-CGA1000	Nonfluorescing longpass filter, transmits >1000 nm	red	green
OF1-CGA760	Nonfluorescing longpass filter, transmits >760 nm	blue	black
OF1-CGA780	Nonfluorescing longpass filter, transmits >780 nm	white	yellow
OF1-CGA830	Nonfluorescing longpass filter, transmits >830 nm	green	orange

4. **Collimating Mirror (specify Standard or SAG+):** The collimating mirror is matched to the 0.22 numerical aperture of our standard optical fibers. Light reflects from this mirror, as a collimated beam, toward the grating. The unit comes with a NIR-enhancing but UV absorbing SAG+ mirror.
5. **Focusing Mirror (specify standard or SAG+):** This mirror focuses first-order spectra on the detector plane. Both the collimating and focusing mirrors are made in-house to guarantee the highest reflectance and the lowest stray light possible. The unit has a standard or SAG+ mirror. As with the collimating mirror, the mirror type needs to be specified when ordering.
6. **Detector:** The detector for Flame-NIR+ is the Hamamatsu G8160-03 InGaAs 128 pixel linear PDA with an effective range of 970-1700nm. The optics split the light into its component wavelengths which fall across the different pixels. Each pixel responds to the wavelength of light that strikes it. The detector outputs an analog signal from each pixel that is converted via the ADC into a digital signal. The driver electronics process this signal and send the spectrum via the USB connection to the software.

Specification	Flame-NIR
Detector	Hamamatsu G8160-03 InGaAs 128 pixel linear PDA
Strengths	Designed as a low cost solution for spectroscopic applications detecting NIR wavelengths
Watch for	Noise levels will increase with elevated temperatures and baseline will rise with long integration times >2 – 3 seconds

See Chapter 6: [Technical Specifications](#) for more detailed detector specifications.

Chapter 6: Technical Specifications

<i>Specification</i>	<i>Flame-NIR</i>
<i>Optical and Spectroscopic</i>	
Integration Time	1 ms to 65 seconds ¹
Dynamic Range for single acquisition ² Typical Observed	7200: 1
Signal-to-Noise (single acquisition)	6000:1
Resolution (FWHM)	10 nm for 25 micron slit
Range	11nm FWHM with 25um slit
Scan rate (max) ⁴	970-1700nm
Spectrometer Channels	One
Thermal Stability	0.08 nm/C (0.014 pixel/C)
Triggering	4 modes (see Triggering Modes)
Triggering Jitter	5 µs
<i>Detector</i>	
Type	Hamamatsu G8160-03 InGaAs linear image sensor
Detector range	900–1700 nm
Pixels	128 pixels
Pixel size	50 µm x 250 µm

<i>Specification</i>	<i>Flame-NIR</i>
Invalid pixels	Minimum:0 Maximum: 2
Electronic shutter	Yes
Pixel well depth	~200,000,000 electrons
Readout noise (single dark spectrum)	9 counts (typical)
Corrected linearity ⁵	>99.8%
Filters (optional)	OF1-CGA830 (preconfigured)
<i>Electrical</i>	
Power requirement (spectrometer functions)	250 mA at +5 VDC
Supply voltage	4.75 – 5.25 V
Power-up time	~2s
Connectors	Micro-USB and JAE DD4 (DD4RA40JA1) 40-pin connector
Micro-USB Absolute Maximum Ratings: V _{CC}	+ 5.5 VDC
DD4 Absolute Maximum Ratings: V _{CC} (Pin 40) Voltage on any pin (other than input power)	+ 5.5 VDC +4VDC
Interface: USB RS-232	USB 2.0, 480 Mbps 2-wire RS-232
<i>Mechanical</i>	
Spectrometer Design	Asymmetric crossed Czerny-Turner
Input Fiber Connector	SMA 905 or FC
Gratings	Grating # NIR33

<i>Specification</i>	<i>Flame-NIR</i>
Entrance Slit	5, 10, 25, 50, 100, or 200 μm slits. (Slits are optional. In the absence of a slit, the fiber acts as the entrance slit.)
Physical Dimensions	88.9 mm x 63.5 mm x 31.9 mm (excluding feet)
Weight	265 g
<i>Environmental</i>	
Temperature: Storage	-30° to +70° C
Operation	0 to 50° C
Humidity	0% - 90% noncondensing
<i>Compliance⁶</i>	
Electrical	CE, FCC, CISPR 11:2010, EMC 2004/108/EC and EN 61326-1:2013
Material	RoHS
Shock	IEC 60068-2-64
Vibration	IEC 60068-2-31
Manufacturing	ISO:9001
¹ Because the baseline will rise rapidly with temperature and extended integration times Ocean Optics recommends integration time be kept below 3 seconds when using the Flame-NIR.	
² Dynamic range for a single acquisition is a measure of the ratio of full signal to noise.	
³ Dynamic range of the system is the range of the detectable light level and can be thought of as the maximum detectable light level at the minimum integration time divided by the minimum detectable light level at the maximum integration time.	
⁴ Scan rate is dependent on the operating computer and not the spectrometer. These figures assume a non real-time operating system.	

<i>Specification</i>	<i>Flame-NIR</i>
⁵ Linearity is measured at 25°C using the linearity correction dark subtraction and the electronic dark correction (if applicable).	
⁶ Contact info@oceanoptics.com to obtain copies of certifications	

Mechanical Diagram


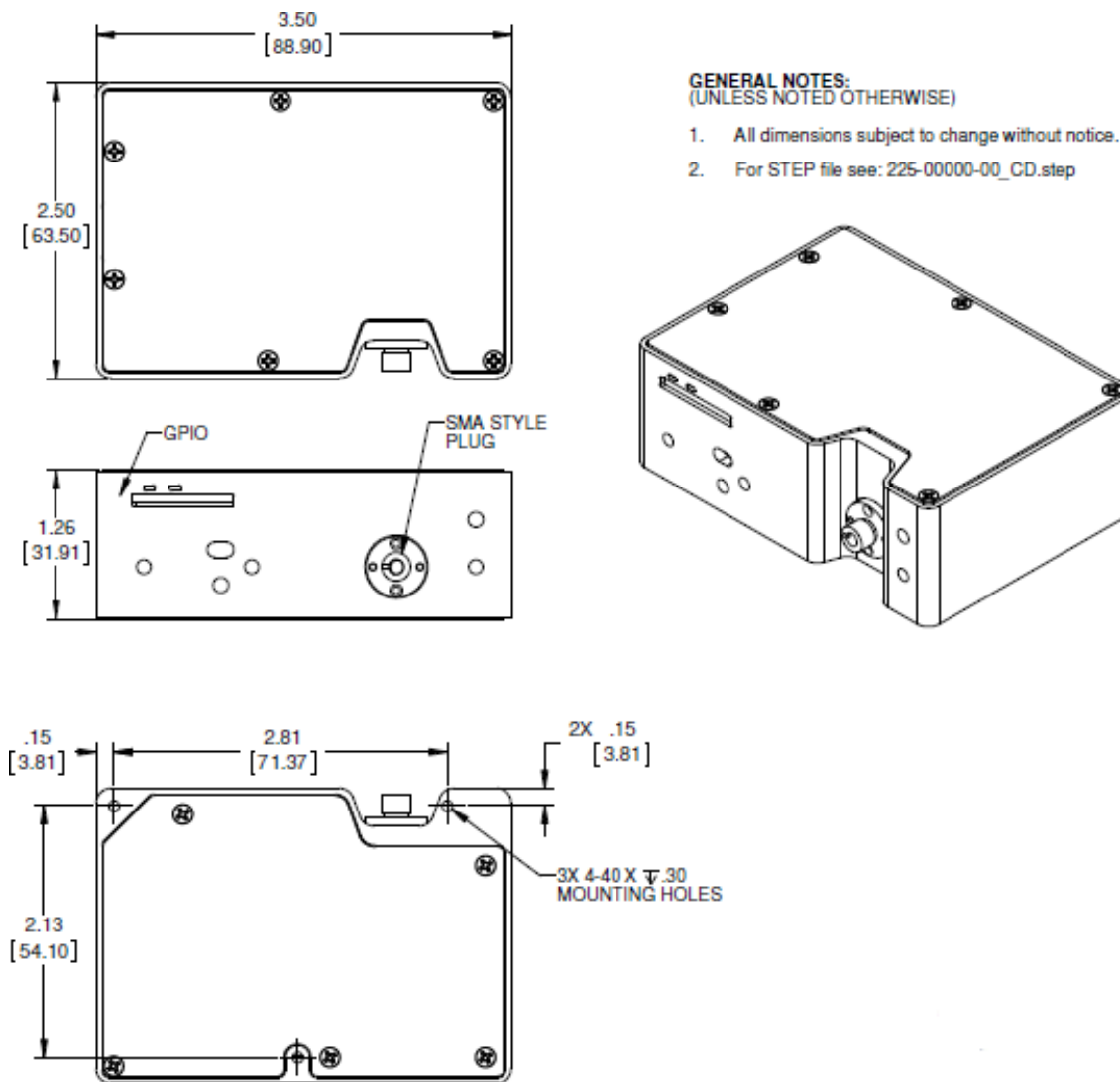


Figure 1: Flame-NIR+ Outer Dimensions

Electrical Pinout

The Flame-NIR+ features a 40-pin Accessory Connector, located on the front of the unit as shown:

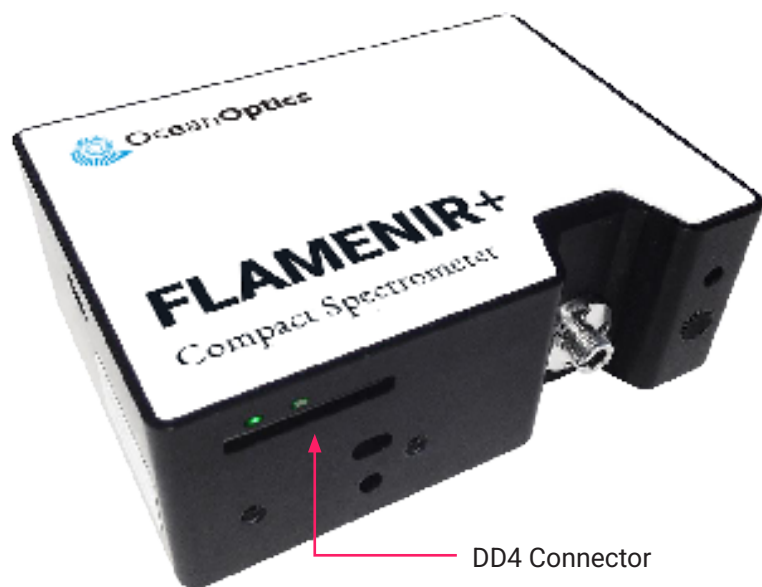

Location of Flame-NIR+ Accessory Connector

DD4 Accessory Connector Pinout Diagram

When facing the 40-pin Accessory Connector on the front of the vertical wall of the Flame-NIR+, pin number 1 is on the right.

Listed below is the pin description for the Flame-NIR+ Accessory Connector located on the front vertical wall of the unit. The Flame-NIR+ will include a JAE DD4 receptacle, part number DD4RA40JA1. Most accessories that plug into the Flame-NIR+ will include a JAE DD4 plug, part number DD4PA40MA1. There is also a vertical connector, JAE part number DD4BA40WA1.

Pin #	Function	Voltage Level	Description
1	Ground	N/A	Ground
2	Trigger	5 or 3.3 V	The TTL trigger signal.
3	Continuous Strobe	5 V	TTL output signal used to pulse a strobe that is divided down from the Master Clock signal.
4	Single Strobe	5 V	TTL output pulse used as a strobe signal, which has a programmable delay relative to the beginning of the spectrometer integration period.
5	Lamp Enable	5 V	A TTL signal that is driven Active HIGH when the Lamp Enable command is sent to the Flame.

6	GPIO 0	2.5V	General Purpose Software Programmable Digital Inputs/Output*
7	GPIO 1	2.5V	General Purpose Software Programmable Digital Inputs/Output*
8	GPIO 2	2.5V	General Purpose Software Programmable Digital Inputs/Output*
9	GPIO 3	2.5V	General Purpose Software Programmable Digital Inputs/Output*
10	Ground	2.5V	General Purpose Software Programmable Digital Inputs/Output*
11	GPIO 4	2.5V	General Purpose Software Programmable Digital Inputs/Output*
12	GPIO 5	2.5V	General Purpose Software Programmable Digital Inputs/Output*
13	GPIO 6	2.5V	General Purpose Software Programmable Digital Inputs/Output*
14	GPIO 7	2.5V	General Purpose Software Programmable Digital Inputs/Output*
15	Ground	N/A	Ground
16	SPI Master Clock	3.3V	Master clock. See SPI below.
17	SPI Master MOSI	3.3V	The SPI Master Out Slave In (MOSI) signal for communications to other SPI peripherals. See SPI below.
18	SPI Master CS	3.3V	TTL output signal used to pulse a strobe that is divided down from the Master Clock signal. See SPI below.
19	SPI Master MISO	3.3V	The SPI Master In Slave Out (MISO) signal for communications to other SPI peripherals. See SPI below.
20	Ground	N/A	Ground
21	I2C Master Clock	3.3V	I2C Master Clock. See I2C below.
22	I2C Master Data	3.3V	I2C Master Data. See I2C below.
23	Ground	N/A	Ground
24	RS232 TX	-6 to +6 V	RS232 Transmit signal for communication with PC connect
25	RS232 RX	N/A	RS232 Receive signal for communication with PC connect
26	Reserved	N/A	Reserved

27	Reserved	N/A	Reserved
28	Reserved	N/A	Do not connect
29	Reserved	N/A	Reserved
30	Reserved	N/A	Reserved
31	Reserved	N/A	Reserved
32	Reserved	N/A	Reserved
33	Reserved	N/A	Reserved
34	Reserved	N/A	Reserved
35	Reserved	N/A	Reserved
36	Reserved	N/A	Reserved
37	Reserved	N/A	Reserved
38	5V Out	5V	The input power pin from the Flame.
39	Ground	N/A	Ground
40	5V In	N/A	The input power pin to the Flame. Additionally when operating via a Universal Serial Bus (USB) this is the USB power connection (+5V) which can be used to power other peripherals (Care must be taken to insure that the peripheral complies with USB Specifications). The entire assembly should not draw more than 500 mA. NOTE: Do not connect both USB power and Auxiliary power (as an input) at the same time.
<i>*See the Caution below.</i>			

CAUTION



Do not connect the GPIO pins to 5V. The GPIOs are not 5V tolerant and will be damaged if connected to 5V. The maximum voltage is 4V for the 3.3V logic pins and 3V for the 2.5V logic pins.

SPI

The Flame-NIR+ has the ability to function as a SPI master through the SPI port, which comprises the SPI Master Clock, SPI Master MOSI, SPI Master CS, and SPI Master MISO pins. To send messages over the SPI port, use the General SPI Input/Output message. The Flame-NIR+ does not send or receive any SPI data without direction from its host PC.



Because SPI is a full-duplex transaction, the General SPI Input/Output message both reads and writes at the same time. For instance, a four byte write will return four bytes of dummy read data, and a four byte read requires four bytes of dummy write data:

- MOSI data is established just prior to the rising edge of the SPI clock
- MISO data is sampled just after a falling edge of the SPI clock.

I2C

The Flame-NIR+ has the ability to function as an I2C master through the I2C port, which comprises the I2C-SDA, and I2C-SCL pins. To send messages over the I2C port, use the General I2C Write and General I2C Read messages. Note that the Flame-NIR+ does not send or receive any I2C data without direction from its host PC. The I2C lines are pulled up internally to 3.3V by 2K resistors.

Performance Charts for Flame-NIR+

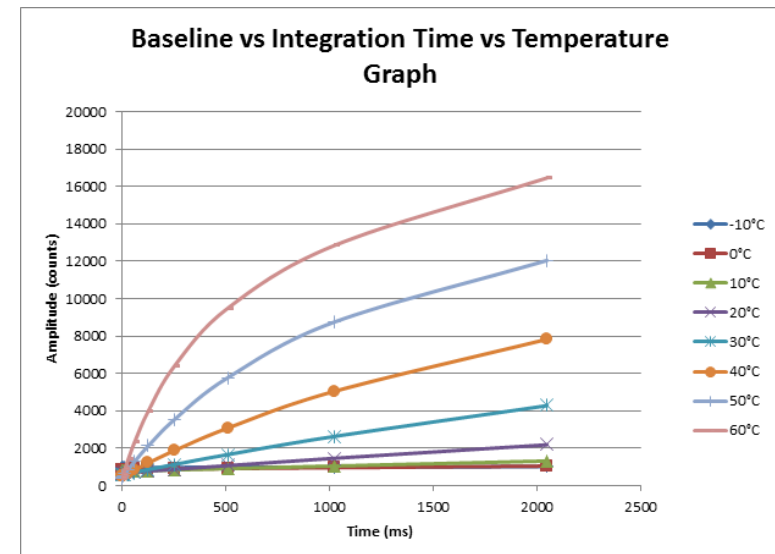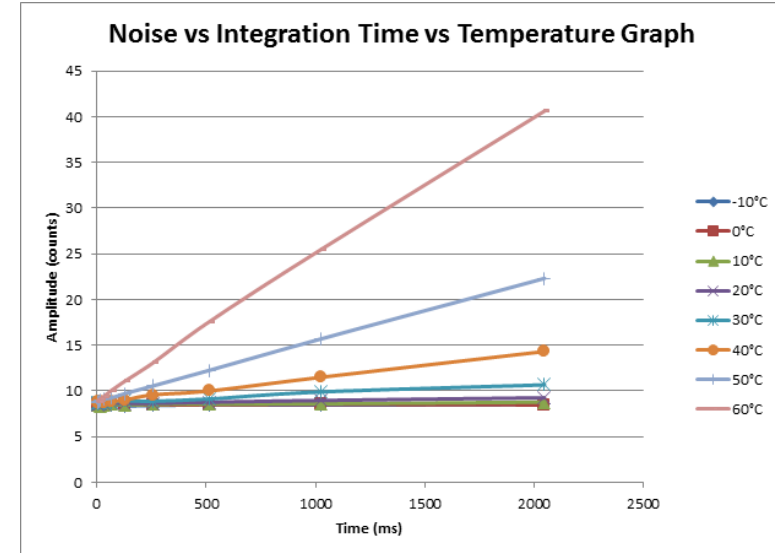

Detector Overview

Detector Description

The detector used for the Flame-NIR+ is an InGaAs photodiode array with integrated CMOS readout circuit.

During each integration period, the current through each photodiode charges an integrated capacitor. After the integration period completes, the charge in each capacitor is measured and then reset. When the capacitor is fully charged, the detector is considered saturated and provides the maximum output level.

Detector Charge Capacity

Ocean Optics products aim for a large signal-to-noise (S:N) ratio (SNR) in optical measurements so that small signal variations can be observed and a large dynamic range is available. The maximum single scan SNR on most CCD based devices is dominated by the well depth. This is because the maximum possible SNR for an integrating photodetector is limited by photon shot noise, which is the square root of the number of electrons collected during the integration period.

The Flame-NIR+, however, has a very large charge storage capacity compared to most CCDs. As a result, the maximum single scan SNR is significantly higher. In fact, photon shot noise is not typically the dominant source of noise in Flame-NIR+ measurements. Instead, SNR is dominated by other sources of noise such as readout noise or thermal noise, depending on the integration time.

Signal Averaging

Signal averaging is an important tool in the measurement of spectral structures. It increases the S:N and the amplitude resolution of a set of samples. The types of signal averaging available in our software are time-based and spatial-based.

When using the time-base type of signal averaging, the S:N increases by the square root of the number of samples. Signal averaging by summing is used

when spectra are fairly stable over the sample period. Thus, even though the maximum single scan SNR is 6000:1 on the Flame-NIR+, a S:N of 60,000:1 should be readily achieved by averaging 100 spectra.

Spatial averaging or pixel boxcar averaging can be used to improve S:N when observed spectral structures are broad. The traditional boxcar algorithm averages n pixel values on each side of a given pixel.

Time-based and spatial-based algorithms are not correlated, so therefore the improvement in S:N is the product of the two processes.

Internal Operation

Pixel Definition

The Flame-NIR+ has 128 pixels, all of which are optically active and not masked. This is slightly different from other Ocean Optics spectrometers which have “dummy” and optical black pixels. Therefore, all pixels read from the Flame NIR in Ocean View, Omnidriver, and directly via USB are active and unmasked. This also means that electronic dark correction is not available on the Flame-NIR+. An invalid pixel is defined as a pixel which has an abnormally high amount of dark current (a “hot” pixel) or which has an abnormally low responsivity (a “dead” pixel)”.

CCD Detector Reset Operation

At the start of each integration period, the detector transfers the signal from each pixel to the readout registers and resets the pixels. The total amount of time required to perform this operation is ~8 - 9 μ s. The user needs to account for this time delay when the pixels are optically inactive, especially in the external triggering modes.

Timing Signals

Strobe Signals

Single Strobe

The Single Strobe signal is a programmable TTL pulse that occurs at a user-determined time during each integration period. This pulse has a user-defined High Transition Delay and Low Transition Delay. The pulse width of the Single Strobe is the difference between these delays. It is only active if the Lamp Enable command is active.

Synchronization of external devices to the spectrometer's integration period is accomplished with this pulse. The Strobe Delay is specified by the Single Strobe High Transition Delay (SSHTD) and the Pulse Width is specified by the Single Strobe Low Transition Delay (SSLTD) minus the Single Strobe High Transition Delay ($PW = SSLTD - SSHTD$). Both values are programmable in 500ns increments for the range of 0 to 65,535 (32.7675ms).

The timing of the Single Strobe is based on the Start of Integration (SOI). SOI occurs on the rising edge of $\overline{X}$ ROG which is used to reset the detector. In all trigger modes using an External Trigger, there is a fixed relationship between the trigger and the SOI. In the Normal mode and Software Trigger mode, the SOI still marks the beginning of the Single Strobe, but due to the nondeterministic timing of the software and computer operating system, this timing will change over time and is not periodic. That is, at a constant integration time, the Single Strobe will not be periodic, but it will indicate the start of the integration. The timing diagram for the Single Strobe in External Hardware Trigger mode is shown below:

The Trigger Delay (TD) is another user programmable delay which specifies the time in 500ns increments that the SOI will be delayed beyond the normal Start of Integration Delay (SOID).

An example calculation of the Single Strobe timing follows:

If the TD = 1ms, SSHTD = 50ms, and SSLTD = 70ms then, the rising edge of the Single Strobe will occur approximately 51.82ms ($1ms + 50ms + 8.2\mu s$) after the External Trigger Input goes high and the Pulse Width will be 20ms ($70ms - 50ms$).

Continuous Strobe

The Continuous Strobe signal is a programmable frequency pulse-train with a 50% duty cycle. It is programmed by specifying the desired period whose range is 2 μ s to 60s. This signal is continuous once enabled, but is not synchronized to the Start of Integration or External Trigger Input. The Continuous Strobe is only active if the Lamp Enable command is active.

Synchronizing Strobe Events

If the application requires more than one pulse per integration period, the user needs to insure the continuous strobe and integration period are synchronized. The integration time must be set so that an equal number of strobe events occurs during any given integration period.

External Triggering

The Flame-NIR+ Spectrometer has several ways of acquiring data. In the Normal/Free-Run mode, the spectrometer is "free running." That is, the spectrometer is continuously scanning, acquiring, and transferring data to your computer, according to parameters set in the software. In this mode, there is no way to synchronize the scanning, acquisition, and transfer of data with an external event. However, trigger pulses for synchronizing an external event with the spectrometer are available.

Each trigger mode involves connecting an external triggering device to the spectrometer and then applying an external trigger to the spectrometer before the software receives the data. The length of the integration time and the source for the integration clock depend upon the mode chosen. All other acquisition parameters are set in the software.


You can trigger the Flame-NIR+ using a variety of External Triggering options through the 40-pin Accessory Connector on the spectrometer.

Triggering Modes

The Flame-NIR+ supports three triggering modes, (plus Normal mode), which are set with the Trigger Mode command.

Normal

In the Normal (Free-run) mode, the spectrometer will acquire one spectrum for each “get spectrum” command issued from the host software. The integration time for each spectrum is pre-programmed prior to the host issuing its “get spectrum” command. The spectrometer waits in a repetitive idling loop until it receives the next “get spectrum” command. Once the command is received, the spectrometer exposes the detector for the pre-programmed integration time, and its resulting spectrum is written to an internal FIFO. The spectrometer signals the host that a spectrum is ready, and the host reads out the new spectrum. Upon reading out the new spectrum, the host immediately commands the spectrometer to acquire the next spectrum. In this process, the spectrometer is responding “on-demand” to the host’s “get spectrum” commands. The spectrometer waits in an idle state between each spectrum and the next “get spectrum” command. The spectrometer does not pre-fetch any spectra in anticipation of the next “get spectrum” command.

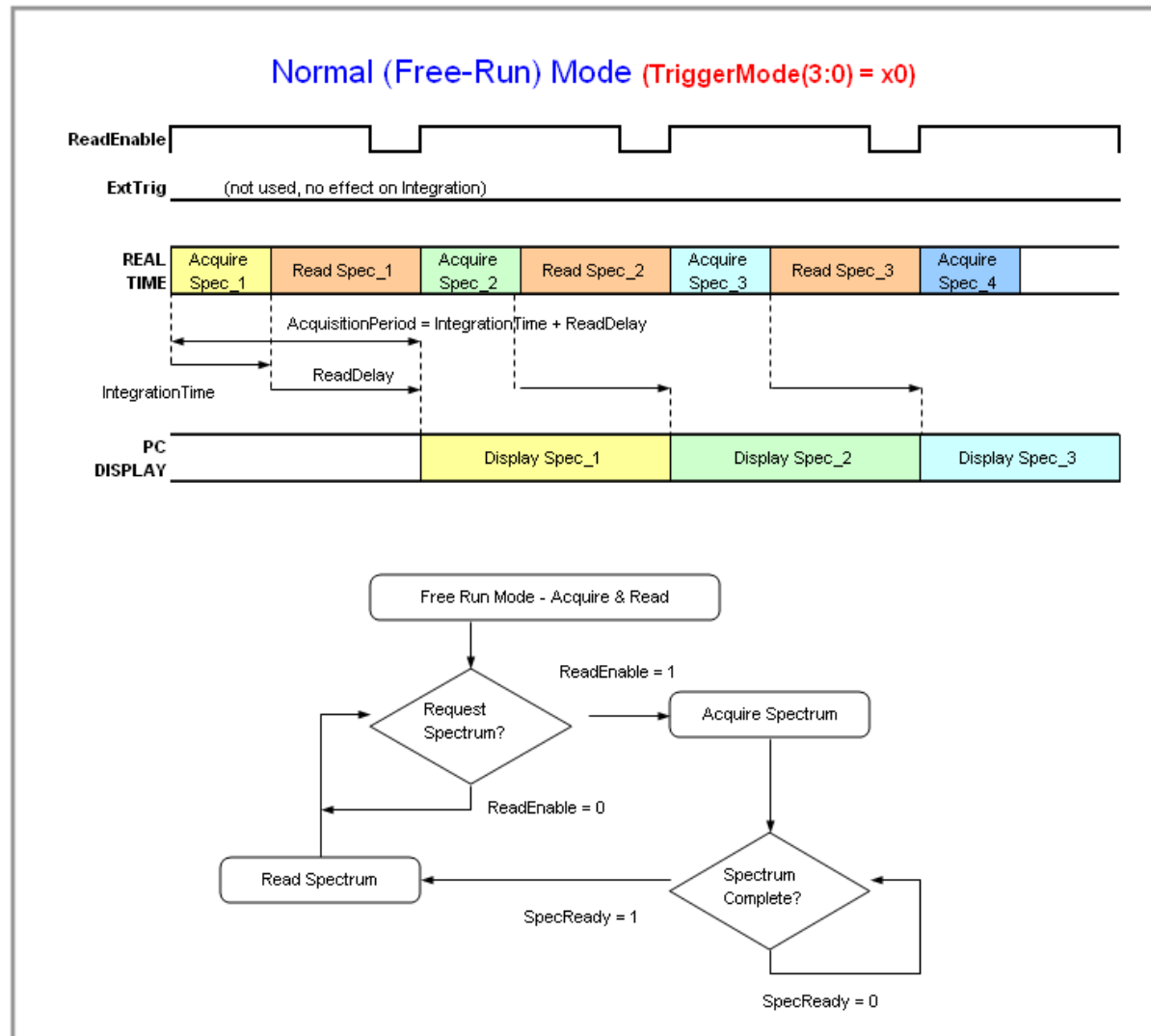


Normal Mode Timing Sequence

External Synchronous Trigger Mode

In the External Synchronous Trigger mode, two external triggers are required to complete a data acquisition. The first rising edge starts the integration period and the second rising edge stops the integration and starts the next. Thus the integration time is the period between the two external trigger pulses. After the integration period, the spectrum is retrieved and available to the user. As in normal mode, no further spectra are acquired until the original spectrum is read by

the user.

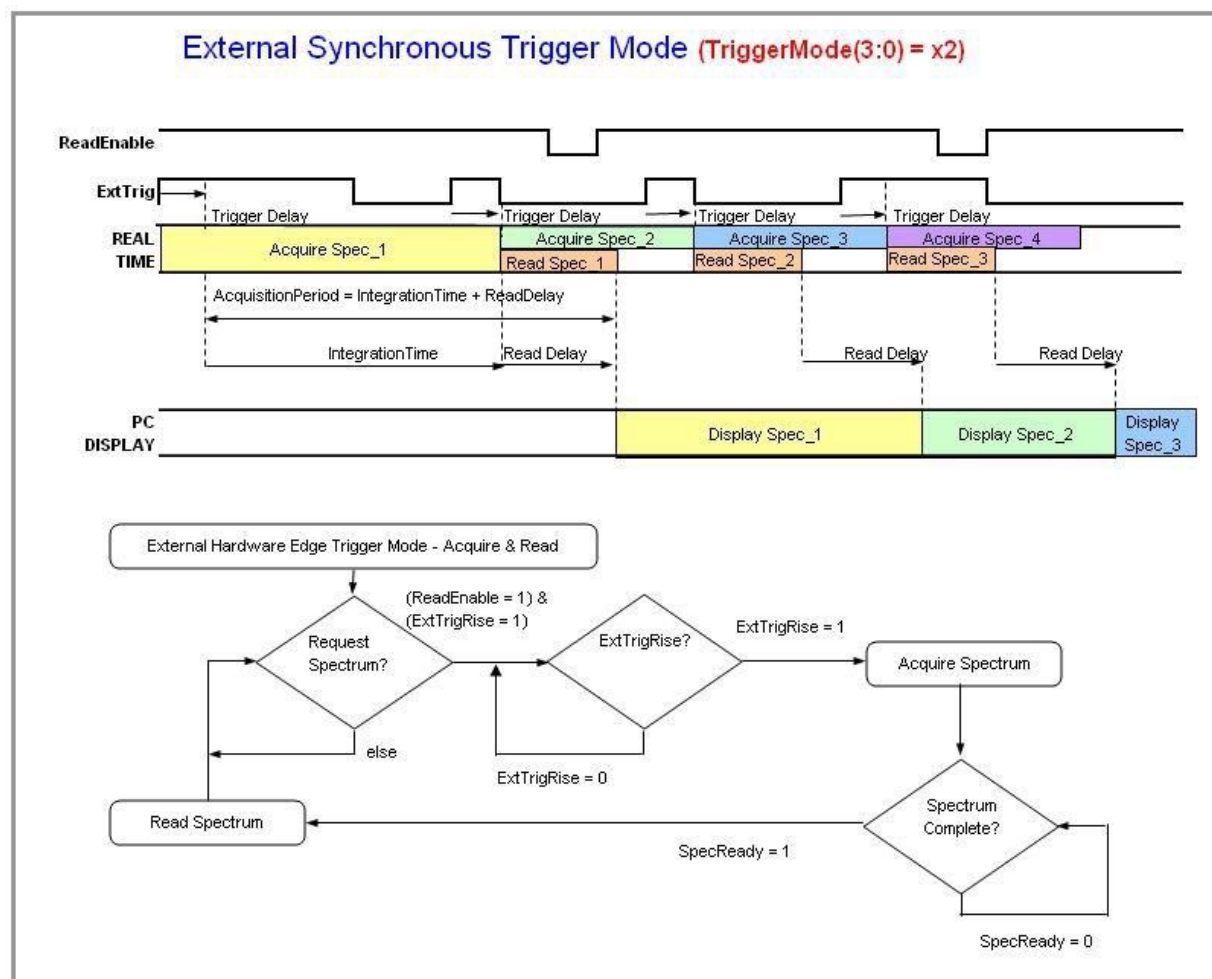


External Synchronous Trigger Mode Timing Sequence

External Hardware Level Trigger Mode

In the External Hardware Level Trigger mode, a rising edge detected by the spectrometer from the External Trigger input starts the integration period specified through the software interface. After the integration period, the spectrum is retrieved and is ready to be read by the user. As long as the trigger level remains active in a logic one state, back-to-back acquisitions can occur, as in the Normal mode, until the trigger transitions to an inactive level. As in normal mode, no

further spectra are acquired until the original spectrum is read by the user.

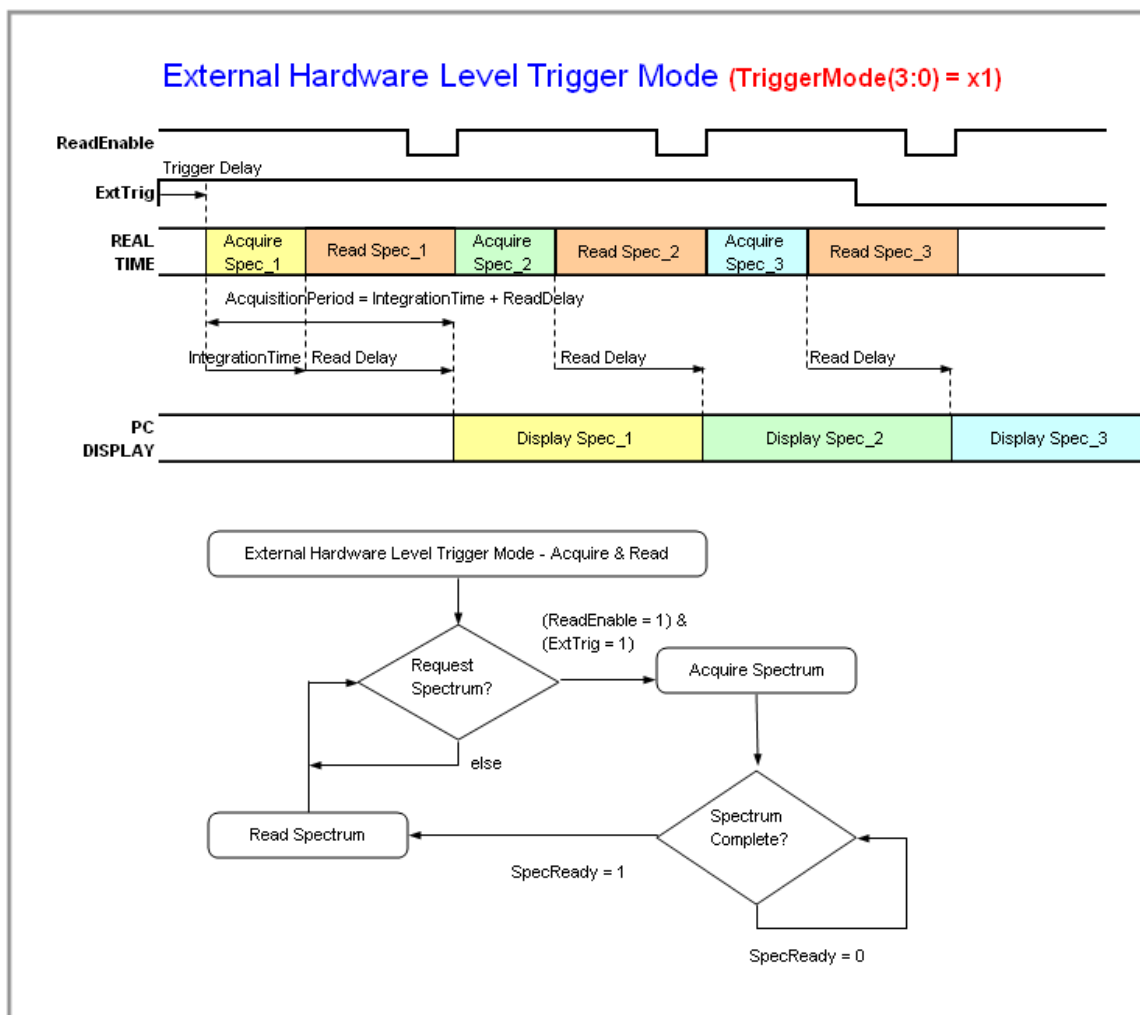


External Hardware Level Trigger Mode Timing Sequence

External Hardware Edge Trigger Mode

In the External Hardware Edge Trigger mode, a rising edge detected by the spectrometer from the External Trigger input starts the integration period specified through the software interface. After the integration period, the spectrum is retrieved and is ready to be read by the user. If another trigger is sent a new integration cycle will begin. If a spectrum request is not received before the integration cycle has ended then that data will be deleted and a new trigger and

spectrum request is required. Only one acquisition will be performed for each External Trigger pulse, no matter what the pulse's duration is. No further spectra are acquired until the original spectrum is read by the user.

HR2000+, USB2000+, Flame-S and -- Timetable for Hardware Edge Trigger Mode
Flame-NIR

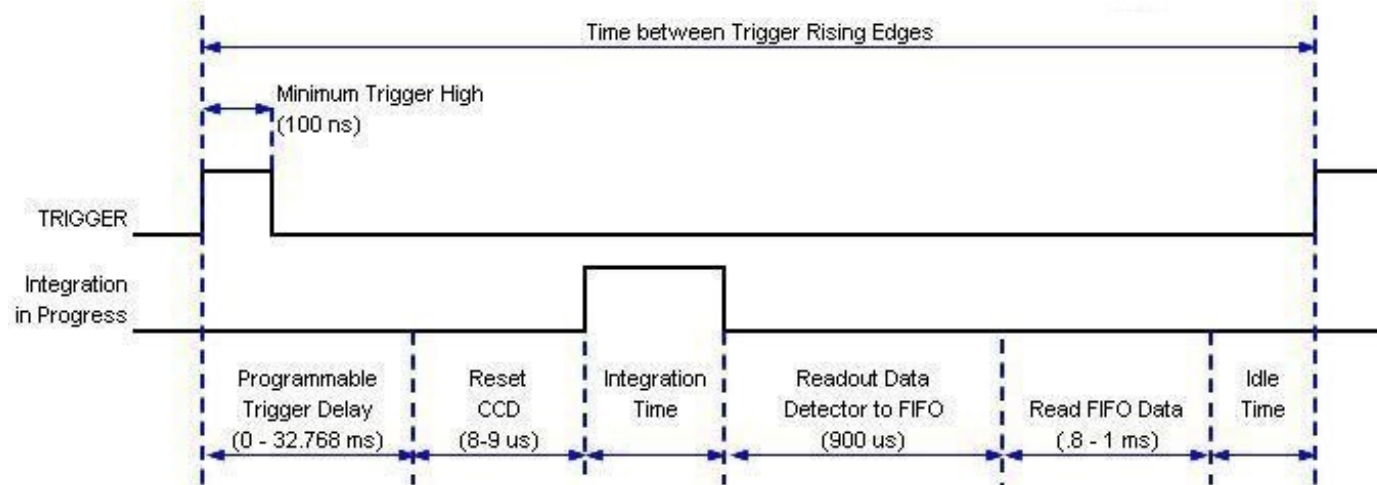

Additional Delay	0 us
Initialize CCD Binning	9 us
Integration time	1,000 us
Read Detector	900 us
Minimum Idle States	1,000 us
Min Trigger Cycle	<u>2,909 us</u>
Min Trigger Cycle	2,909 us

External Hardware Level Trigger Mode Timing Sequence

See DD4 Accessory Connector Pinout Diagram to locate the pins to set up triggering.

Chapter 7: Calibration

Overview

This chapter provides information for performing your own wavelength calibration and irradiance calibration.

An EEPROM flash memory chip in each Flame-NIR+ contains wavelength calibration coefficients, linearity coefficients, and a serial number unique to each individual spectrometer. The spectrometer operating software application reads these values directly from the spectrometer, enabling the ability to “hot-swap” spectrometers between computers without entering the spectrometer coefficients manually on each computer.

Wavelength Calibration

This section describes how to calibrate the wavelength of your spectrometer. Though each spectrometer is calibrated before it leaves Ocean Optics, the wavelength for all spectrometers will drift slightly as a function of time and environmental conditions. Ocean Optics recommends periodically recalibrating the Flame-NIR+.

About Wavelength Calibration

You are going to be solving the following equation, which shows that the relationship between pixel number and wavelength is a third-order polynomial:

$$\lambda_p = I + C_1 p + C_2 p^2 + C_3 p^3$$

Where:

λ = the wavelength of pixel p I = the wavelength of pixel 0

C1 = the first coefficient (nm/pixel)

C2 = the second coefficient (nm/pixel²)

C₃ = the third coefficient (nm/pixel³)

You will be calculating the value for I and the three Cs.

Calibrating the Spectrometer Wavelength

Preparing for Calibration

To recalibrate the wavelength of your spectrometer, you need the following components:

- A light source capable of producing spectral lines

Note: Ocean Optics' HG-2 Mercury-Argon lamp is ideal for recalibration. If you do not have an HG-2, you need a light source that produces several (at least 4-6) spectral lines in the wavelength region of your spectrometer.

- A Flame-NIR+ spectrometer
- An optical fiber (for spectrometers without a built-in slit, a 50-μm fiber works best)
- A spreadsheet program (Excel or Quattro Pro, for example) or a calculator that performs third-order linear regressions

Note: If you are using Microsoft Excel, choose *Tools / Add-Ins* and check *AnalysisToolPak and AnalysisToolPak-VBA*.

Calibrating the Wavelength of the Spectrometer

- **Procedure**

Perform the steps below to calibrate the wavelength of the spectrometer:

1. Place the spectrometer operating software into Quick View (Scope) mode and take a spectrum of your light source. Adjust the integration time (or the A/D conversion frequency) until there are several peaks on the screen that are not off-scale.
2. Move the cursor to one of the peaks and position the cursor so that it is at the point of maximum intensity.
3. Record the pixel number that is displayed in the status bar or legend (located beneath the graph). Repeat this step for all of the peaks in your spectrum.
4. Use the spreadsheet program or calculator to create a table like the one shown in the following figure. In the first column, place the exact or true wavelength of the spectral lines that you used.

In the second column of this worksheet, place the observed pixel number. In the third column, calculate the pixel number squared, and in the fourth column, calculate the pixel number cubed

Independent Variable		Dependent Variables			Values Computed from the Regression	
True Wavelength (nm)		Pixel #	Pixel # ²	Pixel # ³	Predicted Wavelength	Difference
					h	
253.65		175	30625	5359375	253.56	0.09
296.73		296	87616	25934336	296.72	0.01
302.15		312	97344	30371328	302.40	-0.25
313.16		342	116964	40001688	313.02	0.13
334.15		402	161604	64964808	334.19	-0.05
365.02		490	240100	117649000	365.05	-0.04
404.66		604	364816	220348864	404.67	-0.01
407.78		613	375769	230346397	407.78	0.00
435.84		694	481636	334255384	435.65	0.19
546.07		1022	1044484	1067462648	546.13	-0.06
576.96		1116	1245456	1389928896	577.05	-0.09
579.07		1122	1258884	1412467848	579.01	0.06
696.54		1491	2223081	3314613771	696.70	-0.15
706.72		1523	2319529	3532642667	706.62	0.10
727.29		1590	2528100	4019679000	727.24	0.06
738.40		1627	2647129	4306878883	738.53	-0.13
751.47		1669	2785561	4649101309	751.27	0.19

5. Use the spreadsheet or calculator to calculate the wavelength calibration coefficients. In the spreadsheet program, find the functions to perform linear regressions.
 - If using Quattro Pro, look under **Tools / Advanced Math**
 - If using Excel, look under **Analysis ToolPak**
6. Select the true wavelength as the dependent variable (Y). Select the pixel number, pixel number squared, and the pixel number cubed as the independent variables (X). After executing the regression, you will obtain an output similar to the one shown below. Numbers of importance are noted.

Regression Statistics

Multiple R 0.999999831

R Square 0.999999663 ← R Squared

Adjusted R Square 0.999999607

Standard Error 0.125540214

Observations 22

	Intercept	
	Coefficients	Standard Error
Intercept	190.473993	0.369047536
X Variable 1	0.36263983	0.001684745
First coefficient	↑	
X Variable 2	-1.174416E-05	8.35279E-07
Second coefficient	↑	
X Variable 3	-2.523787E-09	2.656608E-10
Third coefficient	↑	

7. Record the Intercept, as well as the First, Second, and Third Coefficients. Additionally, look at the value for R squared. It should be very close to 1. If not, you have most likely assigned one of your wavelengths incorrectly.

Irradiance Calibrations

Irradiance calibrations and relative irradiance calibrations are about quantifying the spectra, by translating the signal (incident number of photons) to a calibration. This can be either absolute (an atomic emission light source of known output power) or relative (corrected for instrument response function but not absolute units). It can be considered a measurement technique and is used widely in remote sensing, light metrology and anywhere where you wish to characterize the incident light source. Irradiance calibrations are not required for many techniques because these measure the relative signal changes with respect to the sample and not the light source.

OceanView has wizards that will step you through absolute irradiance and relative calibrations and more information on these is located in the OceanView manual (see Product-Related Documentation).

Chapter 8: Firmware and Advanced Communications

Flame-NIR+ Firmware

The Flame-NIR+ is a microcontroller-based Miniature Fiber Optic Spectrometer that can communicate via the Universal Serial Bus or RS-232. This section contains the necessary command information for controlling the Flame-NIR+ via the USB interface command. This information is mainly relevant to those who don't wish to use Ocean Optics OmniDriver or SeaBreeze device drivers, or to those who wish to communicate via RS-232.

Hardware Description

The Flame-NIR+ utilizes a Cypress CY7C68013 microcontroller that has a high speed 8051 combined with an USB2.0 ASIC. Program code and data coefficients are stored in external E2PROM that are loaded at boot-up via the I2C bus. The microcontroller has 16K of internal SRAM and 64K of external SRAM. Maximum throughput for spectral data is achieved when data flows directly from the external FIFO's directly across the USB bus. In this mode the 8051 does not have access to the data and thus no manipulation of the data is possible.

USB Information

Ocean Optics Vendor ID number is 0x2457 and the Product ID is 0x104B.

Instruction Set

Command Syntax

The list of the commands is shown in the following table followed by a detailed description of each command. The length of the data depends on the command. All commands are sent to the Flame-NIR+ through End Point 1 Out (EP1). All spectra data is acquired through End Point 2 In and all other queries are retrieved through End Point 1 In (EP1). The endpoints enabled and their order is:

Flame-NIR+ Endpoints

Pipe #	Description	Type	Hi Speed Size (Bytes)	Full Speed Size (Bytes)	Endpoint Address
0	End Point 1 Out	Bulk	512	64	0x01
1	End Point 2 In	Bulk	512	64	0x82
2	End Point 6 In	Bulk	512	64	0x86
3	End Point 1 In	Bulk	512	64	0x81

USB Command Summary

<i>EP1 Command Byte Value</i>	<i>Description</i>
0x01	Initialize Flame-NIR+
0x02	Set Integration Time
0x03	Set Strobe Enable Status
0x04	Set Shutdown Mode
0x05	Query Information
0x06	Write Information

0x09	Request Spectra
0x0A	Set Trigger Mode
0x0B	Query number of Plug-in Accessories Present
0x0C	Query Plug-in Identifiers
0x0D	Detect Plug-ins
0x12	LED Status
0x60	General I ² C Read
0x61	General I ² C Write
0x62	General SPI I/O
0x6A	Write Register Information
0x6B	Read Register Information
0x6C	Read PCB Temperature
0x6D	Read Irradiance Calibration Factors
0x6E	Write Irradiance Calibration Factors
0xFE	Query Information

USB Command Descriptions

A detailed description of all Flame-NIR+ commands follows. While all commands are sent to EP1 over the USB port, the byte sequence is command dependent. The general format is the first byte is the command value and the additional bytes are command specific values.

Byte 0	Byte 1	Byte 2	...	Byte n-1
Command Byte	Command Specific	Command Specific	...	Command Specific

Initialize Flame

Initializes certain parameters on the Flame-NIR+ and sets internal variables based on the USB communication speed the device is operating at. This command should be called at the start of every session; however, if the user does not call it, it will be executed on the first Request Scan command. The default values are set as follows:

Parameter	Default Value
Trigger Mode	0 – Normal Trigger

Byte Format

Byte 0
0x01

Set Integration Time

Sets the Flame-NIR+ integration time in microseconds. The value is a 32-bit value whose acceptable range is 1,000µs – 65,535,000µs. If the value is outside this range the value is unchanged. For integration times less than 655,000µs, the integration counter has a resolution of 10µs. For integration times greater than this the integration counter has a resolution of 1ms.

Byte Format

Byte 0	Byte 1	Byte 2	Byte 3	Byte 4
0x02	LSW-LSB	LSW-MSB	MSW-LSB	MSW-MSB

MSW & LSW: Most Least Significant Word

MSB & LSB: Most/Least Significant Byte

MSW & LSW: Most/Least Significant Word

MSB & LSB: Most/Least Significant Byte

Configuration Index - Description	
0	Serial Number
1	0 th order Wavelength Calibration Coefficient
2	1 st order Wavelength Calibration Coefficient
3	2 nd order Wavelength Calibration Coefficient
4	3 rd order Wavelength Calibration Coefficient
5	Stray light constant
6	0 th order non-linearity correction coefficient
7	0 1 st order non-linearity correction coefficient
8	2 nd order non-linearity correction coefficient
9	3 rd order non-linearity correction coefficient
10	4 th order non-linearity correction coefficient
11	5 th order non-linearity correction coefficient
12	6 th order non-linearity correction coefficient
13	7 th order non-linearity correction coefficient
14	Polynomial order of non-linearity calibration
15	Optical bench configuration: gg fff sss
gg	Grating #, fff – filter wavelength, sss – slit size
16	lame-NIR configuration: AWL V
A	Array coating Mfg, W – Array wavelength (VIS, UV, OFLV), L – L2 lens installed, V – CPLD Version
17	Reserved
18	Reserved
19	29 – User configured

Byte Format

Byte 0	Byte 1
0x05	Configuration Index

Return Format (EP1)

The data is returned in ASCII format and read in by the host through End Point 1.

Byte 0	Byte 1	Byte 2	Byte 3	...	Byte 16
0x05	Configuration Index	ASCII byte 0	ASCII byte 1	...	ASCII byte 14

Write Information

Writes any of the 19 stored spectrometer configuration variables to EEPROM. The 19 configuration variables are indexed as described in the Query Information. The information to be written is transferred as ASCII information.

Byte Format

Byte 0	Byte 1	Byte 2	Byte 3	...	Byte 16
0x06	Configuration Index	ASCII byte 0	ASCII byte 1	...	ASCII byte 14

Request Spectra

Initiates spectra acquisition. The Flame-NIR+ will acquire a complete spectrum. The data is returned in bulk transfer mode through EP2. The table below provides the pixel order overview for the 2 different speeds. The pixel values are decoded as described below.

Byte Format

Byte 0
0x09

Return Format

The format for the returned spectral data is dependent upon the USB communication. The format for both High Speed (480 Mbps) and Full Speed (12Mbps) is shown below. All pixel values are 16 bit values which are organized in LSB | MSB order.

USB High Speed (480Mbps) Packet Format

The data is read from EP2In. The packet format is described below.

Packet #	End Point	# Bytes	Pixels
0	EP2In	256	0-127

The format for the first packet is as follows (all other packets except the synch packet has a similar format except the pixel numbers are incremented by 256 pixels for each packet).

Packet 0

Byte 0	Byte 1	Byte 2	Byte 3
Pixel 0 LSB	Pixel 0 MSB	Pixel 1 LSB	Pixel 1 MSB
...			
		Byte 254	Byte 255
		Pixel 127 LSB	Pixel 127 MSB

Autonulling

Slot 0x11 (17) contains autonulling information that has a scaling term used to adjust the magnitude of the entire spectrum. This can be read out by sending bytes 0x05 11 to the out endpoint (0x01) and then reading out 17 bytes from the in endpoint (0x81). The bytes of use are Byte offset 6 and 7. The 17 bytes will be formatted as follows:

0x05 11 XX XX XX XX SS SS XX XX XX XX XX XX XX XX

Where:

XX = reserved bytes (most are either unused or are only used internally to the device)

SS = saturation level of the device as two bytes (LSB followed by MSB).

These need to be assembled into a single 16-bit value. Any time that a spectrum is read from the spectrometer, each pixel's intensity value should be multiplied by (65535.0/saturation_level) to set the scale appropriately.

The contents of slot 0x11 are set at the factory and should not be altered.

Set Trigger Mode

Sets the Flame-NIR+ Trigger mode to one of four states. If an unacceptable value is passed then the trigger state is unchanged (refer to the External Triggering Options Instructions for Firmware Version 3.0 and Above in Product

Data Value = 0 → Normal (Free running) Mode
Data Value = 1 → External Hardware Level Trigger Mode Data
Value = 2 → External Synchronization Trigger Mode Data
Value = 3 → External Hardware Edge Trigger Mode

Byte Format

Byte 0	Byte 1	Byte 2
0x0A	Data Value LSB	Data Value MSB

Query Number of Plug-in Accessories

To query the number of preset plug-in accessories, use the Query Plug-in Identifiers command 0x0C (below), matching plug-in IDs to the known IDs.

Query Plug-in Identifiers

Queries the Plug-in accessories identifiers. This command returns 7 bytes with the last byte always being zero at this point. Each of the first 6 bytes correspond to Ocean Optics compatible devices which responded appropriately for I2C addresses 2 through 7 respectively. The I2C addresses are reserved for various categories of devices and the value for each category is shown below. I2C addresses 0-1 are reserved for loading program code from EEPROMS

Byte Format

Byte 0
0x0C

Return Format

The data is returned in Binary format and read in by the host through End Point 7.

Byte 0	Byte 1	...	Byte 5	Byte 6
Value @ I ² C address 2	Value @ I ² C address 3	...	Value @ I ² C address 7	0x00

Detect Plug-ins

Reads all of the plug-in accessories that are plugged into the I2C bus. No data values are returned.

Byte Format

Byte 0
0x0D

LED Status

Sets the Flame-NIR+ LEDs as follows.

Byte Format

Byte 0	Byte 1
0x12	Data byte

General I2C Read

Performs a general purpose read on the I2C pins for interfacing to attached peripherals. The time to complete the command is determined by the amount of data transferred and the response time of the peripheral. The I2C bus runs at 400KHz. The maximum number of bytes that can be read is 61.

Command Byte Format

Byte 0	Byte 1	Byte 2
0x60	I ² C Address	Bytes to Read

Return Byte Format

Byte 0	Byte 1	Byte 2	Byte 3	...	Byte N+3
I ² C Results	I ² C Address	Bytes to Read	Data Byte 0	...	Data byte N

<i>I²C Result Value</i>	<i>Description</i>
0	I ² C bus Idle
1	I ² C bus Sending Data
2	I ² C bus Receiving Data
3	I ² C bus Receiving first byte of string
5	I ² C bus in waiting for STOP condition
6	I ² C experienced Bit Error
7	I ² C experience a Not Acknowledge (NAK) Condition
8	I ² C experienced successful transfer
9	I ² C bus timed out

General I2C Write

Performs a general purpose write on the I2C pins for interfacing to attached peripherals. The time to complete the command is determined by the amount of data transferred and the response time of the peripheral. In all I2C communications, the first byte of the transaction consists of a 7 bit address and a read/write bit. The “address” that is passed as the second byte of the I2C write command is this 7 bit address, which will be shifted 1 bit left and appended with the R/W bit to form the first byte of the I2C write transaction. The I2C bus runs at 400KHz. The results codes are described above.

Command Byte Format

<i>Byte 0</i>	<i>Byte 1</i>	<i>Byte 2</i>	<i>Byte 3</i>	<i>...</i>	<i>Byte N+3</i>
0x61	I ² C Address	Bytes to Write	Data Byte 0	...	Data byte N

Return Byte Format

<i>Byte 0</i>
I ² C Results

General SPI Input/Output

Performs a general-purpose write and read on the SPI bus for interfacing to attached peripherals. The time to complete the command is determined by the amount of data transferred and the response time of the peripheral. Wait at least 10 ms after sending a Write command before reading the Return value. The SPI bus runs at ~125KHz Clock. The maximum number of bytes that can be written or read is 61. During this transfer the SPI Chip Select signal is driven to an active LOW TTL level. Data is transmitted out the MOSI (Master Out Slave In) line on the rising edge of the clock signal. Data is also latched in the from the MISO line on the falling edge of the clock signal.

Command Byte Format

<i>Byte 0</i>	<i>Byte 1</i>	<i>Byte 2</i>	<i>Byte 3</i>	<i>...</i>	<i>Byte N+2</i>
0x62	# of Bytes (N)	Write Byte 0	Write Byte 1	...	Write Byte N

Return Byte Format

<i>Byte 0</i>	<i>Byte 1</i>	<i>Byte 2</i>	<i>...</i>	<i>Byte N+1</i>
# of Bytes (N)	Read Byte 0	Read Byte 1	...	Read Byte N

Write Register Information

Most all of the controllable parameters for the Flame-NIR+ are accessible through this command (e.g., GPIO, strobe parameters, etc.). A complete list of these parameters with the associate register information is shown in the table below. Commands are written to End Point 1 Out typically with 4 bytes (some commands may require more data bytes). All data values are 16 bit

values transferred in LSB | MSB order. This command requires 100µs to complete; the calling program needs to delay for this length of time before issuing another command. In some instances, other commands will also write to these registers (i.e., integration time), in these cases the user has the options of setting the parameters through 2 different methods.

Byte Format

Byte 0	Byte 1	Byte 2	Byte 3
0x6A	Register Value	Data Byte LSB	Data Byte MSB

Register Address	Description	Default Value	Min Value	Max Value	Time Base
0x00*	Master Clock Counter Divisor	24	1	0xFFFF	48MHz
0x04	FPGA Firmware Version (Read Only)				
0x08	Continuous Strobe Timer Interval Divisor	48000	0	0xFFFF	Continuous Strobe Base Clock (see Register 0x0C)
0x0C	Continuous Strobe Base Clock Divisor	4800	0	0xFFFF	48MHz
0x0C	Continuous Strobe LSB Register	4800	0	0xFFFF	48MHz
0x10*	Integration Period Base Clock Divisor	480	0	0xFFFF	48MHz
0x10*	Integration Period LSB Register	480	0	0xFFFF	1KHz
0x14	Set base_clk or base_clkx2 0: base_clk 1: base_clkx2	0	0	1	N/A
0x18*	Integration Clock Timer Divisor	600	0	0xFFFF	Integration Period Base Clock (see Register 0x10)
0x18*	Integration Period MSB Register				
0x20	Reserved				

0x28	Hardware Trigger Delay – Number of Master Clock cycles to delay when in External Hardware Trigger mode before the start of the integration period	0	0	0xFFFF	
0x28	Hardware Trigger Delay – Delay the start of integration from the rising edge of the trigger in 500ns increments	0	0	0xFFFF	

Register Address	Description	Default Value	Min Value	Max Value	Time Base
0x2C*	Trigger Mode 0 = Free Running 1 = External Hardware Level Trigger 1 = External Synchronization Trigger 3 = External Hardware Edge Trigger	0	0	4	N/A
0x30	Reserved				
0x38	Single Strobe High Clock Transition Delay Count	1	0	0xFFFF	2MHz
0x3C	Single Strobe Low Clock Transition Delay Count	5	0	0xFFFF	2MHz
0x40	Lamp Enable	0	0	1	N/A
0x44	High Gain	0	0	1	N/A
0x48	GPIO Mux Register 0: pin is GPIO pin 1: pin is alternate function	0	0	0x03FF	N/A
0x50	GPIO Output Enable 1: pin is output 0: pin is input	0	0	0x03FF	N/A
0x54	GPIO Data Register For Output: Write value of signal For Input: Read current GPIO state	0	0	0x03FF	N/A
0x58	Reserved				

0x5C	Offset Value	0	0	0xFFFF	N/A
0x60	Offset Control Bit 0 = Enable Auto-Nulling Bit 1 = Enable Auto-Nulling Saturation	0	0	0xFFFF	N/A
0x64	FPGA Programmed (Read Only)	0x5501	N/A	N/A	N/A
0x68	Maximum Saturation Level	0x55F0	0	0xFFFF	N/A

Notes:

* - User should not change these values because spectrometer performance can be affected. This information is included just for completeness
& - These values are controlled by other command interfaces to the Flame-NIR+ (i.e., Set integration time command).

Read Register Information

Read the values from any of the registers above. This command is sent to End Point 1 Out and the data is retrieved through End Point 1 In.

Byte Format

Byte 0	Byte 1
0x6B	Register Value

Return Format (EP1In)

Byte 0	Byte 1	Byte 2
Register Value	Value LSB	Value MSB

Read PCB Temperature

Read the Printed Circuit Board Temperature. The Flame-NIR+ contains a DS1721 temperature sensor chip which is mounted to the underside of the PCB. This command is sent to End Point 1 Out and the data is retrieved through End Point 1 In. The value returned is a signed 16-bit A/D conversion value, which is equated to temperature by:

$$\text{Temperature (}^{\circ}\text{C)} = .003906 * \text{ADC Value}$$

Byte Format

Byte 0
0x6C

Return Format (EP1In)

Byte 0	Byte 1	Byte 2
Read Result	ADC Value LSB	ADC Value MSB

If the operation was successful, the Read Result byte value will be 0x08. All other values indicate the operation was unsuccessful.

Read Irradiance Factors

Reads 60 bytes of data, which is utilized for Irradiance Calibration information from the desired EEPROM memory address.

Byte Format

Byte 0	Byte 1	Byte 2
0x6D	EEPROM Address LSB	EEPROM Address MSB

Return Byte Format

Byte 0	Byte 1	...	Byte 59
Byte 0	Byte 1	...	Byte 59

Write Irradiance Factors

Write 60 bytes of data, which is utilized for Irradiance Calibration information to the desired EEPROM memory address.

Byte Format

Byte 0	Byte 1	Byte 2	Byte 3	...	Byte 62
0x6E	EEPROM Address LSB	EEPROM Address MSB	Byte 0	...	Byte 59

Query Status

Returns a packet of information containing the current operating information. Packet structure is given below:

Byte Format

Byte 0
0xFE

Return Format

The data is returned in Binary format and read in by the host through End Point 1 In. The structure for the return information is as follows:

Byte	Description	Comments
0-1	Number of Pixels - WORD	LSB MSB order
2-5	Integration Time - WORD	Integration time in μ s – LSW MSW. Within each word order is LSB MSB

Byte	Description	Comments
6	Lamp Enable	0 – Signal LOW 1 – Signal HIGH
7	Trigger Mode Value	
8	Spectral Acquisition Status	
9	Packets In Spectra	Returns the number of Packets in a Request Spectra Command.
10	Power Down Flag	0 – Circuit is powered down 1 – Circuit is powered up
11	Packet Count	Number of packets loaded into End Point Memory
12	Reserved	
13	Reserved	
14	USB Communications Speed	0 – Full Speed (12Mbps) 0x80 – High Speed (480 Mbps)
15	Reserved	

Flame-NIR+ Serial Port Interface

Communications and Control Information

The Flame-NIR+ is a microcontroller-based Miniature Fiber Optic, which can communicate via the Universal Serial Bus or RS-232. This document contains the necessary command information for controlling the Flame-NIR+ via the RS-232 interface.

Hardware Description

The Flame-NIR+ utilizes a Cypress FX2 microcontroller, which has a high speed 8051, combined with an USB ASIC. Program code and data coefficients are stored in external E2PROM, which are loaded at boot-up via the I2C bus.

Instruction Set

Command Syntax

The list of the commands is shown in the following table along with the microcode version number they were introduced with. All commands consist of an ASCII character passed over the serial port, followed by some data. The length of the data depends on the command. The format for the data is either ASCII or binary (default). The ASCII mode is set with the “a” command and the binary mode with the “b” command. To insure accurate communications, all commands respond with an ACK (ASCII 6) for an acceptable command or a NAK (ASCII 21) for an unacceptable command (i.e., data value specified out of range).

In the ASCII data value mode, the Flame-NIR+ “echoes” the command back out the RS-232 port. In binary mode all data, except where noted, passes as 16-bit unsigned integers (WORDS) with

the MSB followed by the LSB. By issuing the “v command” (Version number query), the data mode can be determined by viewing the response (ASCII or binary).

In a typical data acquisition session, the user sends commands to implement the desired spectral acquisition parameters (integration time, etc.). Then the user sends commands to acquire spectra (S command) with the previously set parameters. If necessary, the baud rate can be changed at the beginning of this sequence to speed up the data transmission process.

RS232 Command Summary

Letter	Description
A	Adds scans
B	Set Pixel Boxcar
C	
D	
E	
F	***Non functional but follows SAD500 command format***
G	Set Data Compression
H	***Non functional but follows SAD500 command format***
I	Sets integration time
J	Sets Lamp Enable Line
K	Changes baud rate
L	Clear Memory
M	Set Data Storage Mode
N	
O	
P	Partial Pixel Mode

Q	
R	
S	Starts spectral acquisition with previously set parameters
T	Sets trigger mode
U	

Letter	Description
V	
W	Query scans in memory
X	
Y	
Z	Read out Scan from memory
A	Set ASCII mode for data values
b	Set binary mode for data values
k	Sets Checksum mode
o	Oxygen (USB-LS-450) related commands
t	Temperature in hundredths of degrees C
u	Set Oxygen Calibration Coefficients
v	Provides microcode version #
x	Sets calibration coefficients
y	Sets 16-bit timer operation

?	Queries parameter values
+	Reads the plugged-in accessories
-	USB2000 Identifier

Command Descriptions

A detailed description of all Flame-NIR+ commands follows. The {} indicates a data value which is interpreted as either ASCII or binary (default). The default value indicates the value of the parameter upon power up.

Add Scans

Sets the number of discrete spectra to be summed together. Since the Flame-NIR+ has the ability to return 32 bit values, overflow of the raw 16 bit ADC value is not a concern.

Command Syntax:	A{DATA WORD}
Response:	ACK or NAK

Range:	1-5000
Default value:	1

Pixel Boxcar Width

Sets the number of pixels to be averaged together. A value of n specifies the averaging of n pixels to the right and n pixels to the left. This routine uses 32-bit integers so that intermediate overflow will not occur; however, the result is truncated to a 16-bit integer prior to transmission of the data. This math is performed just prior to each pixel value being transmitted out. Values greater than ~3 will exceed the idle time between values and slow down the overall transfer process.

Command Syntax: <i>B{DATA WORD}</i>	
Response:	ACK or NAK
Range:	0-15
Default value:	0

Set Data Compression

Specifies whether the data transmitted from the Flame-NIR+ should be compressed to speed data transfer rates.

Command Syntax: <i>G{DATA WORD}</i>	
Response:	ACK or NAK
Range:	0 – Compression off !0 – Compression on
Default value:	0

Integration Time (16 bit)

Sets the Flame's integration time, in milliseconds, to the value specified.

Command Syntax: <i>I{16 bit DATA WORD}</i>	
Response:	ACK or NAK
Range:	1 - 65000
Default value:	10

Integration Time (32 bit)

Sets the Flame's integration time, in microseconds, to the value specified.

Command Syntax: <i>i{32 bit DATA WORD}</i>	
Response:	ACK or NAK
Range:	1000 – 65,000,000
Default value:	10,000

Lamp Enable

Sets the Flame's Lamp Enable line to the value specified

Command Syntax: <i>J{DATA WORD}</i>	
Value:	0 = Light source/strobe off—Lamp Enable low 1 = Light source/strobe on—Lamp Enable high
Response:	ACK or NAK
Default value:	0

Baud Rate

Sets the Flame's baud rate.

Command Syntax: <i>K{DATA WORD}</i>	
Value:	0=2400 1=4800 2=9600 3=19200 4=38400 5=Not Supported 6=115,200
Response:	See below
Default value:	2

When changing baud rates, the following sequence must be followed:

1. Controlling program sends K with desired baud rate, communicating at the old baud rate
2. The Flame-NIR+ responds with ACK at old baud rate, otherwise it responds with NAK and the process is aborted
3. Controlling program waits longer than 50 milliseconds
4. Controlling program sends K, followed by the desired baud rate. For example, to change the baud rate from 9600 to 115,200 in ASCII mode:
 - At 9600 baud, send "K6<enter>"
 - You will receive an ACK at 9600 baud
 - Send "K6<enter>" at 115200. (Note that the "K" is not echoed back, but the "6" is)
 - You will receive two ACK characters and the prompt at 115200 baud.
4. The Flame-NIR+ responds with ACK at new baud rate, otherwise it responds with NAK and old baud rate is used



If a deviation occurs at any step, the previous baud rate is utilized.

Clear Memory

Clears the spectral data memory based on the valued specified. Clearing memory is immediate since only pointer values are reinitialized.

CAUTION

All stored spectra are lost when the Clear memory command is executed.

Command Syntax:	L{DATA WORD}
Value:	0= Clear spectral memory 1= Clear spectral memory
Response:	ACK or NAK
Default value:	N/A

Data Storage Mode

Sets the data storage mode for future spectral acquisitions.

Command Syntax:	M{DATA WORD}
Value:	0= Scans transmitted through the serial port 1= Scans stored in spectral memory and not transmitted
Response:	ACK or NAK
Default value:	0

Pixel Mode

Specifies which pixels are transmitted. While all pixels are acquired on every scan, this parameter determines which pixels will be transmitted out the serial port.

Command Syntax:	P{DATA WORD}	
Value:	Description 0 = all 127 pixels 1 = every nth pixel with no averaging 2 = N/A 3 = pixel x through y every n pixels 4 = up to 10 selected pixels between 0 and 127 (denoted p1, p2, ... p10) 5 = up to 10 pixel ranges between 0 and 2047 (denoted r1start, r1end, r2start, r2end, ...r10start, r10end)	Example P 0 (spaces for clarity only) P 1<Enter> N<Enter> P 2 N/A P3<Enter> x<Enter> y<Enter> n<Enter> P 4<Enter> n<Enter> p1<Enter> p 2<Enter> p3<Enter> ... p10<Enter> P5<Enter> n<Enter> r1start<Enter> r1end<Enter> r2start<Enter> r2end<Enter>...?
Response:	ACK or NAK	
Default value:	0	



Since most applications only require a subset of the spectrum, this mode can greatly reduce the amount of time required to transmit a spectrum while still providing all of the desired data. This mode is helpful when interfacing to PLCs or other processing equipment.

Spectral Acquisition

Acquires spectra with the current set of operating parameters. When executed, this command determines the amount of memory required. If sufficient memory does not exist, an ETX (ASCII 3) is immediately returned and no spectra are acquired. An STX (ASCII 2) is sent once the data is acquired and stored. If the Data Storage Mode value is 0, then the data is transmitted immediately.

Command Syntax:	S
Response:	If successful, STX followed by data If unsuccessful, ETX

The format of returned spectra includes a header and an end of spectrum pattern, in addition to 128 active pixels. The format is as follows:

WORD 0xFFFF – start of spectrum
WORD Data size flag (0=Data is WORD's, 1=Data is DWORD's) WORD Number of Scans Accumulated
WORD Integration time in milliseconds
WORD FPGA Established Baseline value (LSW) WORD FPGA Established Baseline value (MSW) WORD pixel mode
WORDS if pixel mode not 0, indicates parameters passed to the Pixel Mode command (P) (D)WORDS spectral data depending on Data size flag (128 pixels)
WORD 0xFFFD – end of spectrum

Trigger Mode

Sets the Flame's external trigger mode to the value specified.

Command Syntax:	T{DATA WORD}
Value:	0 = Normal (Free running) Mode 2 = External Hardware Level Trigger Mode 3 = External Synchronization Trigger Mode 4 = External Hardware Edge Trigger Mode Note: This is not consistent with the USB trigger modes.
Response:	ACK or NAK
Default value:	0

Set FPGA Register Value

Sets the appropriate register within the FPGA. The list of register setting is in the USB command set information. This command requires two data values, one to specify the register and the next to specify the value.

Command Syntax:	W{DATA WORD 1}{DATA WORD 2}
Value:	Data Word 1 – FPGA Register address Data Word 2 – FPGA Register Value
Response:	ACK or NAK
Default value:	N/A

ASCII Data Mode

Sets the mode in which data values are interpreted to be ASCII. Only unsigned integer values (0 – 65535) are allowed in this mode and the data values are terminated with a carriage return (ASCII 13) or linefeed (ASCII 10). In this mode the Flame-NIR+ "echoes" the command and data values back out the RS-232 port.

Command Syntax:	aA
Response:	ACK or NAK
Default value	N/A



The command requires that the string “aA” be sent without any CR or LF. This is an attempt to insure that this mode is not entered inadvertently.

A legible response to the Version number query (v command) indicates the Flame-NIR+ is in the ASCII data mode.

Binary Data Mode

Sets the mode in which data values are interpreted to be binary. Only 16 bit unsigned integer values (0 – 65535) are allowed in this mode with the MSB followed by the LSB

Command Syntax:	bB
Response:	ACK or NAK
Default value	Default at power up – not changed by Q command



The command requires that the string “bB” be sent without any CR or LF. This is an attempt to insure that this mode is not entered inadvertently.

Checksum Mode

Specifies whether the Flame-NIR+ will generate and transmit a 16-bit checksum of the spectral data. This checksum can be used to test the validity of the spectral data, and its use is recommended when reliable data scans are required.

Command Syntax:	k{DATA WORD}
Value:	0 = Do not transmit checksum value !0 = transmit checksum value at end of scan
Response:	ACK or NAK
Default value:	0

Version Number Query

Returns the version number of the code running on the microcontroller. A returned value of 1000 is interpreted as 1.00.0.

Command Syntax:	v
Response:	ACK followed by {DATA WORD}
Default value	N/A

Calibration Constants

Writes one of the 16 possible calibration constant to EEPROM. The calibration constant is specified by the first DATA WORD which follows the x. The calibration constant is stored as an ASCII string with a max length of 15 characters. The string does not check to see if it makes sense.

Command Syntax:	x{DATA WORD}{ASCII STRING}
Value:	DATA WORD Index description 0 – Serial Number 1 – 0 th order Wavelength Calibration Coefficient 2 – 1 st order Wavelength Calibration Coefficient 3 – 2 nd order Wavelength Calibration Coefficient 4 – 3 rd order Wavelength Calibration Coefficient 5 – Stray light constant 6 – 0 th order non-linearity correction coefficient 7 – 1 st order non-linearity correction coefficient 8 – 2 nd order non-linearity correction coefficient 9 – 3 rd order non-linearity correction coefficient 10 – 4 th order non-linearity correction coefficient 11 – 5 th order non-linearity correction coefficient 12 – 6 th order non-linearity correction coefficient 13 – 7 th order non-linearity correction coefficient 14 – Polynomial order of non-linearity calibration 15 – Optical bench configuration: gg fff sss gg – Grating #, fff – filter wavelength, sss – slit size 16 – Flame-NIR+ configuration: AWL V A – Array coating Mfg, W – Array wavelength (VIS, UV, OFLV), L – L2 lens installed, V – CPLD Version 17 – Reserved 18 – Reserved 19 – Reserved
Response:	ACK or NAK
Default value:	N/A

To query the constants, use the ?x{DATA WORD} format to specify the desired constant

Query Variable

Returns the current value of the parameter specified. The syntax of this command requires two ASCII characters. The second ASCII character corresponds to the command character which sets the parameter of interest (acceptable values are B, A, I, K, T, J, i, j, y). Note that 'i' returns the integration time rounded to the nearest 10us, where 'j' returns 1μsec resolution. A special case of this command is ?x (lower case) which requires an additional data word be passed to indicate which calibration constant is to be queried.

Command Syntax:	?{ASCII character}
Response:	ACK followed by {DATA WORD}
Default value:	N/A

Questions?

Chat with us at www.oceanoptics.com.
info@oceanoptics.com

US +1 727-733-2447

EUROPE +31 26-3190500

ASIA +86 21-6295-6600

We Measure What Matters.

We empower scientists, engineers, and innovators with precise, reliable optical sensing solutions that drive discovery and improve lives.